

## 1 Purpose

This document provides release notes for TID21, a data management software for Metrohm handheld Raman spectrometers including products of TacticID Plus, TacticID 1064, TacticID Mobile and TacticID-1064ST.

The new features, improvements, fixed bugs and known issues for each version of TID21 are included.

For installation refer to **809538101EN TID21 Installation Guide and User Manual** for details.

Release Notes:

- [TID21 v2.01.1.45](#)
- [TID21 v2.01.1.36](#)
- [TID21 v2.01.1.35](#)
- [TID21 v2.01.1.33](#)

## 2 TID21 v2.1.1.45

Details the changes in version 2.01.1.45 over the previously officially released version of 2.01.1.35 or 2.01.1.36.

Release Date: 2025.08



## 2.3 Improvements

- Default database server is named Plainsboro (for new installations).
- Default server instance name changed to *Plainsboro*
- Report preview removed. Export button now immediately launches the export screen.
- Blue and gray added as background colors when editing library samples.
- Update does not require the user to reconfigure the database or restart the server.
- RNDIS driver included in the TID21 installation folder.

## 2.4 Fixed Bugs

- Fixed bug causing device to appear in TID21 on the data transfer screen after manually disconnecting the USB cable.
- Fixed bug causing device's factory number to display instead of the device's serial number.
- Fixed bug causing the Export button to be active while nothing was selected in the Library screen.
- Fixed bug where TID21 was not displaying the user set UDF name in Summary Report.
- Fixed bug where server instance names are not listed under IT.

## 2.5 Known Issues

- n/a.

### 3 TID21 v2.1.1.36

Details the changes in version 2.01.1.36 over the previously officially released version of 2.01.1.27 or 2.01.1.35.

Release Date: 2025.04



## 4.1 Compatibility

To ensure compatibility for software operation, you must use matching versions of TID21 on the PC and TOS on the device.

*Table 4 TID21 v2.1.1.35 Compatibility*

|                              |                  |                 |                     |                                |
|------------------------------|------------------|-----------------|---------------------|--------------------------------|
| <b>Product</b>               | TacticID-1064 ST | TacticID Mobile | TacticID Plus       | TacticID-1064<br>(BWS492-1064) |
| <b>Embedded<br/>firmware</b> | TOS-XM 2.1.2080  | TOS-XM 2.1.1373 | TOS Plus 2.01.227.0 | TOS1064 2.01.227.0             |

It's highly recommended to uninstall old TID software to ensure the data integrity.

## 4.2 New Features

- Added Czech language support.

### 4.3 Improvements

- Login interface for Spanish language fixed.
- Translation errors fixed.

## 4.4 Fixed Bugs

- Fixed bug stopping Users, Verification, and Results screens from being scrolled.

## 4.5 Known Issues

TID21 may crash while accessing records.

## 5 TID21 v2.1.1.33

Details the changes in version 2.01.1.33 over the previously officially released version of 1.01.78.1.

Release Date: 2024.10

