

1 Purpose

This document describes new functions, the changes, improvements and fixed bugs and problems in the following firmware version:

- 57.1007.0012 for OS version 005
- in comparison with the firmware version:
- 57.1007.0011 for OS version 004

2 New functions

The following new functions have been integrated:

- Support for PostScript USB printers
- Support for Sartorius balances via COM port (USB/RS-232 Converter required)
- COM port settings for balances

3 Improvements

The following improvements were made:

- Firmware optimizations for the Russian language.
- Installation wizard is activated at initial instrument start-up.

4 Fixed bugs and problems

The following bugs and problems were fixed:

- Some translation mistakes were corrected.
- OHMDOS-759 – The dosing rate was not updated correctly in the Live Display.

5 Known bugs and problems

The following bugs or problems are known:

- none