



## Application Note AN-EC-037

# Differential electrochemical mass spectrometry

## Integrating Metrohm Autolab potentiostats with Hiden mass spectrometers


Differential electrochemical mass spectrometry (DEMS) is an advanced analytical technique that integrates an electrochemical experiment with mass spectrometry. This method enables real-time, *in situ* analysis of gaseous and volatile electrochemical reactants, intermediates, and products. DEMS is particularly valuable for studying reaction mechanisms and kinetics in energy conversion and

storage applications, such as fuel cells, batteries, and electrocatalytic processes.

This Application Note highlights both the DEMS technique itself and the combination of a Hiden mass spectrometer with a VIONIC used to detect and quantify hydrogen production during a cyclic voltammogram (CV).



Automatic current ranging should also be disabled in order to ensure that the current passed to the Hiden software is correct. Using a trigger cable, INTELLO is able to trigger the mass spectrometer to begin scanning. The Hiden software utilizes a falling edge trigger. The initial/end state of the DIO port on the potentiostat (Figure 3) prior to the measurement should therefore be set to high (1). Pin 5 is used for this.

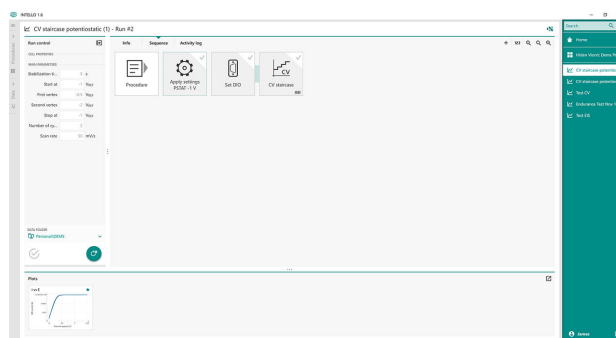

**Figure 3.** End-state of the instrument in INTELLO.

## EXPERIMENTAL


The bespoke cable used in these experiments is shown in Figure 4. The 25-pin side connects to the HPR-40 system. The 15-pin side is for the VIONIC, and there are two BNC connectors for the *I* and *E* signals, respectively.



**Figure 4.** DIO trigger and BNC cable used in this system.

At the beginning of the procedure, the state of pin 5 should be changed to low (0). This is done with a Set DIO command (Figure 5).


All Hiden software has the ability to accept triggers either as a Start/Stop automation event (MASsoft) or in the triggering window (QGA 2). When the trigger signal is received the currently selected setup file will start automatically.

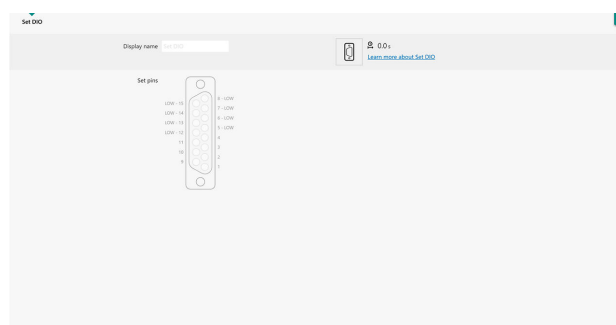

**Figure 5.** Set DIO command detailed view.

## RESULTS AND DISCUSSION


Cyclic voltammetry (Figure 6) was conducted in the region of -0.5 to -2 V, with a start/stop voltage of -1 V vs AgCl. The scan rate was 50 mV/s. There are no discernable features in the CV until about -1.3 V, where the production of hydrogen is triggered via the reduction of the aqueous electrolyte.

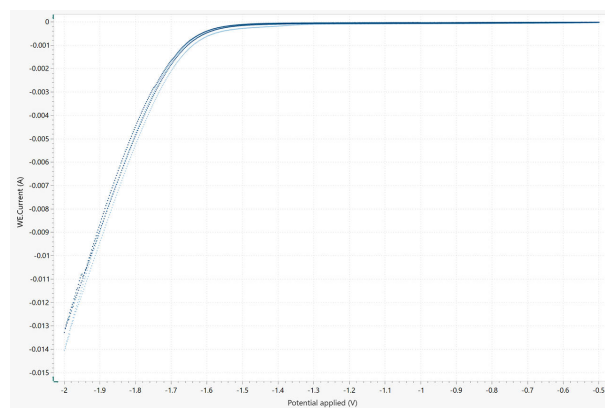

**Figure 6.** Cyclic voltammogram of 1 mol/L NaOH on an Au-sputtered electrode.

The synchronized mass, current, and potential vs time signals are shown in Figure 7. No discernable change in the mass signal is observed until -1.3 V is reached.

The rise in signal corresponds with (and also confirms) the production of hydrogen at the very negative potentials.

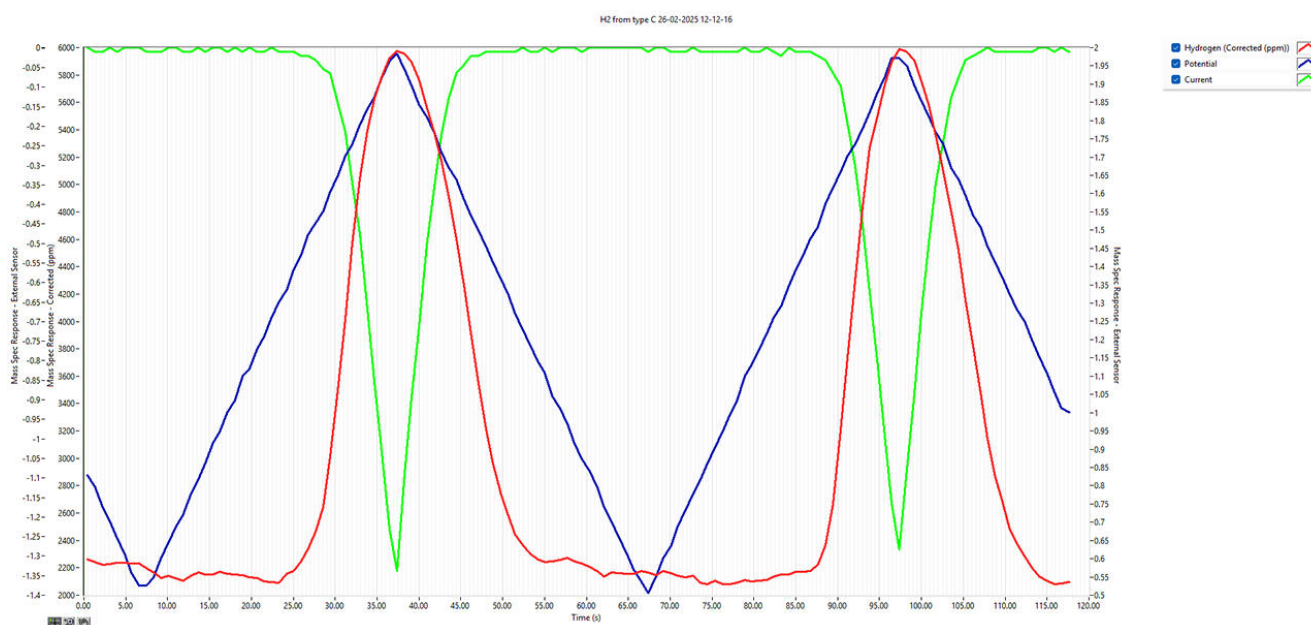

**Figure 7.** Synchronized, mass (red), potential (blue), and current (green) signals from the Hiden software.

## REFERENCE

1. Clark, E. L.; Bell, A. T. Direct Observation of the Local Reaction Environment during the Electrochemical Reduction of CO<sub>2</sub>. *J. Am. Chem. Soc.* **2018**, *140* (22), 7012–7020. <https://doi.org/10.1021/jacs.8b04058>.

## CONTACT

Metrohm Nordic ApS  
Lyskær 3C. 1. th.  
2730 Herlev

[mail@metrohm.dk](mailto:mail@metrohm.dk)

## CONFIGURATION



### VIONIC

VIONIC is our new-generation potentiostat/galvanostat that is powered by Autolab's new INTELLO software.

VIONIC offers the **most versatile combined specifications of any single instrument** currently on the market.

- Compliance voltage:  $\pm 50$  V
- Standard current  $\pm 6$  A
- EIS frequency: up to 10 MHz
- Sampling interval: down to 1  $\mu$ s

Also included in VIONIC's price are features that would usually carry an additional cost with most other instruments such as:

- Electrochemical Impedance Spectroscopy (EIS)
- Selectable Floating
- Second Sense (S2)
- Analog Scan