



Application Note AN-P-086

Quality assurance of instant coffee

Free and total carbohydrate determination with IC-PAD according to AOAC 996.04 and ISO 11292

Coffee is an extremely popular beverage with a significant economic importance. Quality assurance, including tracing adulterants in coffee, is therefore an established process and a requirement for consumers.

Carbohydrates which make up to 50% of raw coffee beans function as flavor, viscosity, and aroma agents [1]. They also serve as authenticity tracers, because unadulterated soluble coffee is exclusively made from pure roasted coffee [2,3]. There are clear specification criteria for the quality assessment by ISO 24114 and AFCASOLE (e.g., a limit of <2.46% total glucose and <0.45%

total xylose expressed as mass fractions of total carbohydrates) [3]. AOAC 996.04 and ISO 11292 give analytical requirements for instant coffee quality tests regarding analysis of free and total carbohydrates. Ion chromatography (IC) allows precise quantification of the mandatory analytes arabinose, fructose, galactose, glucose, mannose, sucrose, mannitol, and xylose according to AOAC and ISO. The presented IC method is extremely sensitive and overcomes a very common challenge of possible analyte coelutions, such as with rhamnose.

SAMPLE AND SAMPLE PREPARATION

Instant coffee powders (300 mg per 100 mL) of two instant coffee brands (Jacobs coffee GOLD and a customer sample) were prepared as described in AOAC and ISO to determine the amount of free carbohydrates (arabinose, fructose, galactose, glucose, mannose, sucrose, and mannitol) and total carbohydrates (arabinose, galactose, glucose, mannose, xylose, and mannitol) after acid hydrolyzation.

For the determination of free carbohydrates, coffee powders were dissolved in 100 mL ultrapure water (UPW) and then filtered (0.25 μm). For total carbohydrate analysis, coffee powders were hydrolyzed in HCl (0.1 mol/L) at 100 ° C (150 minutes), diluted to 100 mL with UPW, and filtered with an $\text{Ag}^+\text{-H}^+$ -cartridge combination. A final dilution (10 to 50-fold) is recommended with UPW.

EXPERIMENTAL

The carbohydrates specified above for dissolved (free) and total carbohydrate analysis were baseline separated on a Metrosep Carb 2 column with a binary high-pressure gradient combined with a flow gradient (940 Professional

IC Vario ONE/HPG configuration) (**Figure 1**). Amperometric detection was performed after PCR with 300 mmol/L NaOH to improve the detection sensitivity of the method.

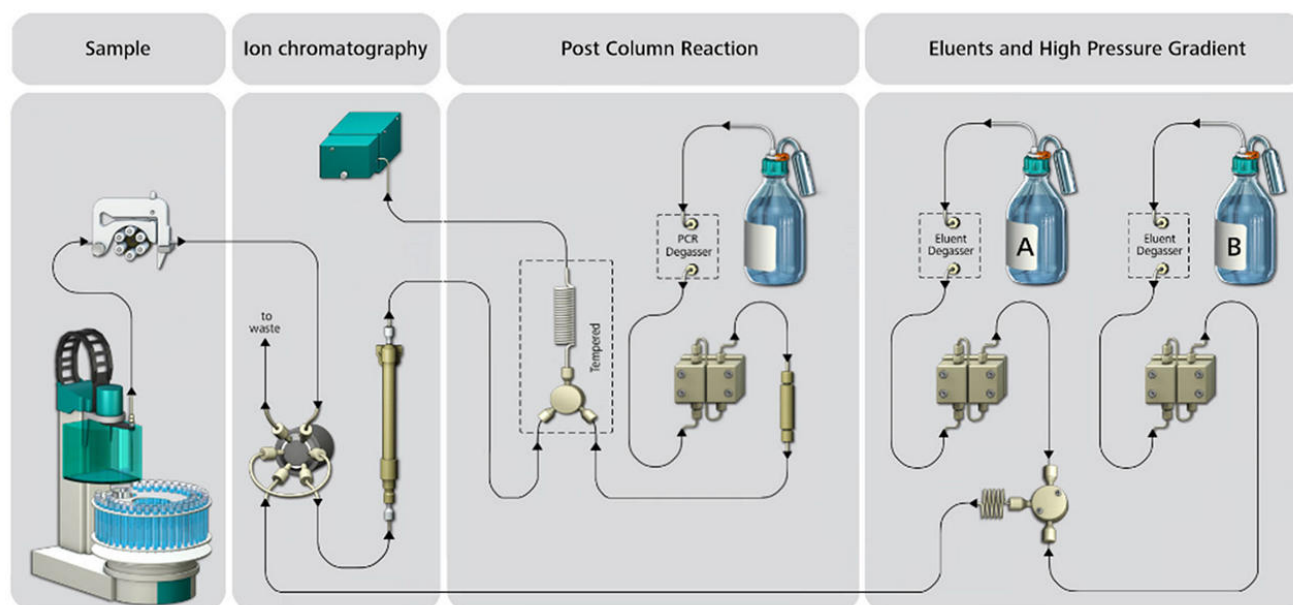


Figure 1. Schematic showing the sample flow path from sample introduction with an 858 Professional Sample Processor, to the 940 Professional IC with column (Metrosep Carb 2 - 250/4.0), amperometric detector (Wall-Jet cell with Au and Pd electrodes), and the high pressure gradient pumps for eluent A (UPW) and B (200 mmol/L NaOH and 1 mmol/L NaAc). To increase sensitivity, 300 mmol/L NaOH is added as a PCR (Post Column Reaction) solution. Chromatography for anions is often referred as HPAEC (high performance anion exchange chromatography) but is simplified here to the generic term of IC.

RESULTS

For the two analyzed instant coffee samples, the free carbohydrate content (results not shown) after dissolution in UPW ranged from 0.2 to 27 g/kg. The mass fractions show unique patterns for both samples. In the Jacobs brand, arabinose and mannose dominate (35 mass%), while the largest peaks for the customer-provided instant coffee brand corresponded to glucose (20 mass%) and fructose (almost 40 mass%).

Total carbohydrate content after acid hydrolysis is especially crucial for quality control and purity assessment (Table 1). ISO 24114 set limitations

for total glucose and xylose of 2.32 and 0.42%, respectively. The total carbohydrate content of both tested samples show a distinct distribution (Table 1 and Figure 2). Both brands contain similar fractions of galactose and arabinose. Glucose, mannose, and xylose contents differ in a broader range.

Looking more closely at the quality criteria, the purity of Jacobs coffee GOLD can be approved as unadulterated product. The customer-provided brand indicates adulteration and would fail a respective control.

Table 1. Carbohydrate concentrations (g/kg) determined by IC-PCR- PAD after acid hydrolysis in two instant coffee samples (Jacobs coffee GOLD and a customer sample). The total carbohydrate content is expressed as the individual mass fractions (M%) of mannitol, arabinose, galactose, glucose, mannose, and xylose (ISO 11292). Additionally, quantification of rhamnose, fructose, ribose and sucrose is possible (Figure 2). Purity indicators are given by the limits for total glucose (<2.32%) and total xylose (<0.42%) (ISO 24114:2011).

	Jacobs (g/kg) [M%]	Customer (g/kg) [M%]
Mannitol	ND	9 [2.6%]
Arabinose	28.3 [6.5%]	36 [10.2%]
Galactose	190.0 [43.9%]	197.8 [56.2%]
Glucose	6.3 [1.5%]	34.2 [9.7%]
Mannose	207.1 [47.8%]	68.5 [19.4%]
Xylose	1.2 [0.3%]	6.7 [1.9%]
Total carbohydrate content	436.9 [100%]	352.2 [100%]

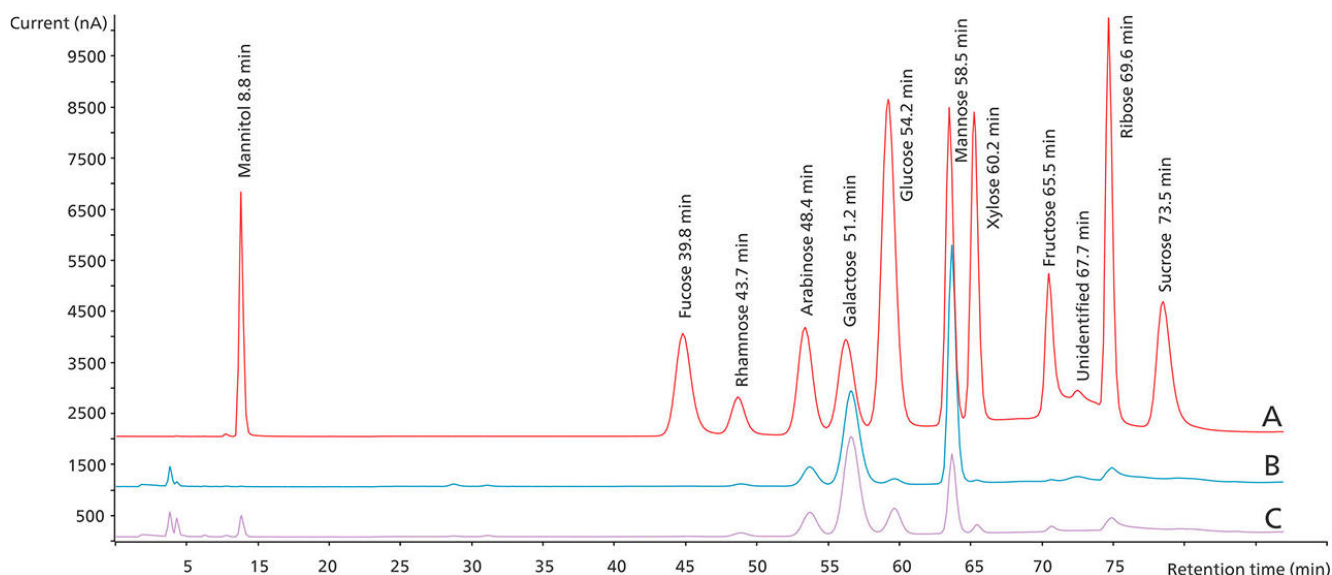


Figure 2. Chromatogram overlay of (A) a 5 mg/kg mixed carbohydrate standard and the diluted (1:10) samples of (B) Jacobs coffee and (C) customer instant coffee after acid hydrolysis. The separation and detection were performed according to the setup listed in Figure 1. For better performance, quantification of fructose and ribose should be performed using peak height.

CONCLUSION

With the presented method the requirements of **AOAC 996.04** and **ISO 11292** for the determination of dissolved and total carbohydrates in instant coffee are fulfilled. An excellent separation of the required carbohydrates can be achieved by combination of a binary high-pressure gradient and a flow gradient on a Metrosep Carb 2 column. An additional benefit of this method eliminates peak overlap between rhamnose and arabinose, an overall constraint of the ISO-method. Overall,

the precise quantification of all required carbohydrates plus fucose and ribose can be performed.

Automation and Inline Sample Preparation are additional improvements to increase the sample throughput and save laboratory time and money.

IC with amperometric detection is a robust, highly specific and precise valuable addition for analytical laboratories performing carbohydrate analysis.

REFERENCES

1. Araya and Rao (2007), Crit Rev Food Sci Nutr. 47(1), 51–67.
2. Girard et al. (2006), J AOAC Int. 89(4), 99–1003.
3. AFCASOLE (Association of European producers of soluble coffee) statement on the authenticity of soluble coffees of 6 July 1995; as confirmed by the ECF (European Coffee Federation, legal successor of AFCASOLE) in January 2007.

CONTACT

メトロームジャパン株式会
社
143-0006 東京都大田区平
和島6-1-1
null 東京流通センター アネ
ックス9階

metrohm.jp@metrohm.jp

CONFIGURATION




940 Professional IC Vario ONE/HPG

940 プロフェッショナル IC Vario ONE/HPG はバイナリー高圧グラジエントを備えたサプレッションの無しのインテリジェントコンパクトIC装置です。942 拡張モジュール を使用することでクォーターグラジエントシステムにまで拡張することができます。この装置は任意の分離メソッドおよび検出メソッドによって使用することができます。

典型的な使用領域:

- グラジエント溶離法後のパルスタイプのアンペロメトリック検出器による炭水化物分析
- ポストカラム法を使用する、もしくは使用しない UV/VIS検出器によるグラジエント使用

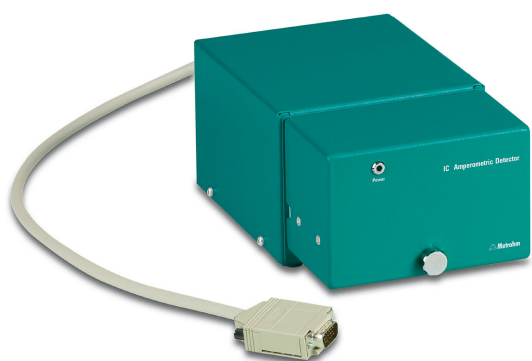

IC Amperometric Detector

インテリジェントIC装置のためのコンパクトかつインテリジェントなアンペロメトリック検出器。4つの測定モード(DC、PAD、flexIPAD、CV)による優れた選択性、卓越した信号・騒音比、非常に高速の測定待機機能が高精度の測定できます。



Metrosep Carb 2 - 250/4.0

ICカラムMetrosep Carb 2 - 250/4.0は、アルカリ性溶離液およびパルスアンペロメトリー検出を用いた炭水化物の測定に特に適しています。大容量の陰イオン交換カラムは、スチロール/ジビニルベンゼン共重合体をベースとしています。これはpH = 0~14の範囲で安定しており、単糖類および二糖類を分離します。さらに、数種の糖アルコール、無水糖、アミノ糖などの分析にも適しています。Metrosep Carb 2分離カラムの250mmのバージョンは、複雑な分離のために最適化されています。



858 Professional Sample Processor – Pump

858 プロフェッショナルサンプルプロセッサ - ポンプは、500 μ Lから500 mLまでのサンプルを処理します。サンプルは内蔵式双方向性の2チャンネルのペリスタリックポンプまたは800 ドジーノ電動ビュレットによって転送されます。



942 Extension Module Vario HPG

942 拡張モジュール Vario HPGは、プロフェッショナル IC Vario装置群のための拡張モジュールです。このモジュールにより、プロフェッショナルIC Vario高圧グラジエントシステムで更にもう1種類の溶離液使用することが可能になります。

典型的な使用領域:

- 最大4種類の溶離液までの高圧グラジエントシステム



IC equipment Wall-Jet cell: Carb (Au, Pd)

Equipment comprised of Wall-Jet cell with additional accessories. For carbohydrate analysis with a gold working electrode and a Pd reference electrode.