Application Note AN-S-404

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1. U.S. EPA. U.S. EPA Method 300.1:
Determination of Inorganic Anions in
Drinking Water by Ion Chromatography,
Revision 1.0, 1997.
2. Boorman, G. A.; Dellarco, V.; Dunnick, J. K.;
et al. Drinking Water Disinfection
Byproducts: Review and Approach to
Toxicity Evaluation. *Environmental Health
Perspectives* **1999**, *107*, 207–217.
[DOI:10.2307/3434484](https://doi.org/10.2307/3434484)
3. Jackson, P. E. Ion Chromatography in
Environmental Analysis. In *Encyclopedia of
Analytical Chemistry*; Meyers, R. A., Ed.;
Wiley, 2000.
[DOI:10.1002/9780470027318.a0835](https://doi.org/10.1002/9780470027318.a0835)
4. ISO. ISO 10304-4:2022 - Water Quality —
Determination of Dissolved Anions by
Liquid Chromatography of Ions — Part 4:
Determination of Chlorate, Chloride and
Chlorite in Water with Low
Contamination, 2022.
5. ISO. ISO 10304-1:2007 - Water Quality —
Determination of Dissolved Anions by
Liquid Chromatography of Ions — Part 1:
Determination of Bromide, Chloride,
Fluoride, Nitrate, Nitrite, Phosphate and
Sulfate, 2007.

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Metrosep A Supp 21 - 250/4.0

The Metrosep A Supp 21 columns are designed for operation with hydroxide-based eluents and provide excellent separating efficiency, coupled with a very high capacity. The small particles (4.6 μm) based on hydrophilized polystyrene/divinylbenzene guarantee sharp peaks. The stationary phase exhibits high stability with respect to temperature, pressure, and pH value, and is therefore suitable for extreme working conditions. The longer Metrosep A Supp 21 - 250/4.0 column version was specially developed for the determination of oxohalides (chlorite, bromate, chlorate), standard anions (fluoride, chloride, nitrite, bromide, nitrate, sulfate, and phosphate), and DCAA (dichloroacetate). With its separating efficiency, it exceeds the requirements of the US EPA method 300.1 A+B and of the DIN EN ISO 10304-1&4 standard. The high column capacity enables the quantification of anions and oxohalides in low $\mu\text{g/L}$ concentrations with excellent reproducibility, even in the most challenging sample matrices. With the wide range of elution conditions available, it is also possible to determine other anionic components, e.g. low-molecular-weight organic acids.

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