Application Note AN-COR-016

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REFERENCES

1. G61 Standard Test Method for Conducting Cyclic Potentiodynamic Polarization Measurements for Localized Corrosion Susceptibility of Iron-, Nickel-, or Cobalt-Based Alloys.
<https://www.astm.org/standards/g61>
(accessed 2024-05-24).
2. Standard Practice for Conventions Applicable to Electrochemical Measurements in Corrosion Testing.
<https://www.astm.org/g0003-14r19.html>
(accessed 2024-03-08).
3. Bellezze, T.; Viceré, A.; Giuliani, G.; et al. Study of Localized Corrosion of AISI 430 and AISI 304 Batches Having Different Roughness. *Metals* **2018**, 8 (4), 244.
[DOI:10.3390/met8040244](https://doi.org/10.3390/met8040244)

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metrohm.jp@metrohm.jp

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Platinum sheet electrode

This electrode consists of a thin sheet of platinum embedded into a glass shaft. This electrode can be used as a counter electrode for most electrochemical measurements. The surface area is approximatively 1 cm².

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1 L corrosion cell

The Autolab 1 L corrosion cell is suitable for corrosion measurements according to ASTM standards. The cell has a thermostatic jacket for temperature control and a series of openings for counter electrodes, pH sensor, thermometer, Luggin-Haber capillary and gas purging. The 1 l corrosion cell has been designed to measure the corrosion properties samples with a diameter ranging from 14.7 mm to 16 mm and a thickness from 0.5 mm up to 4 mm. The exposed surface is 1 cm² and the seals are made out of natural rubber.

VIONIC

VIONIC is our new-generation potentiostat/galvanostat that is powered by Autolab's new INTELLO software. VIONIC offers the most versatile combined specifications of any single instrument currently on the market. Compliance voltage: ± 50 V; Standard current ± 6 A; EIS frequency: up to 10 MHz; Sampling interval: down to 1 μ s; Also included in VIONIC's price are features that would usually carry an additional cost with most other instruments such as: Electrochemical Impedance Spectroscopy (EIS); Selectable Floating; Second Sense (S2); Analog Scan;