Application Note AN-EC-033

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1. Morávková, Z.; Dmitrieva, E. Structural Changes in Polyaniline near the Middle Oxidation Peak Studied by in Situ Raman Spectroelectrochemistry. *Journal of Raman Spectroscopy* **2017**, 48 (9), 1229–1234. <https://doi.org/10.1002/jrs.5197>.
2. Dong, J.-C.; Zhang, X.-G.; Briega-Martos, V.; et al. In Situ Raman Spectroscopic Evidence for Oxygen Reduction Reaction Intermediates at Platinum Single-Crystal Surfaces. *Nat Energy* **2019**, 4 (1), 60–67. <https://doi.org/10.1038/s41560-018-0292-z>.
3. Lopez-Ramirez, M. R.; Aranda Ruiz, D.; Avila Ferrer, F. J.; et al. Analysis of the Potential Dependent Surface-Enhanced Raman Scattering of p-Aminothiophenol on the Basis of MS-CASPT2 Calculations. *J. Phys. Chem. C* **2016**, 120 (34), 19322–19328. <https://doi.org/10.1021/acs.jpcc.6b05891>.
4. Tabatabaei, M.; Sangar, A.; Kazemi-Zanjani, N.; et al. Optical Properties of Silver and Gold Tetrahedral Nanopyramid Arrays Prepared by Nanosphere Lithography. *J. Phys. Chem. C* **2013**, 117 (28), 14778–14786. <https://doi.org/10.1021/jp405125c>.

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metrohm.jp@metrohm.jp

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VIONIC

VIONIC is our new-generation potentiostat/galvanostat that is powered by Autolab's new INTELLO software. VIONIC offers the most versatile combined specifications of any single instrument currently on the market. Compliance voltage: ± 50 V; Standard current ± 6 A; EIS frequency: up to 10 MHz; Sampling interval: down to 1 μ s; Also included in VIONIC's price are features that would usually carry an additional cost with most other instruments such as: Electrochemical Impedance Spectroscopy (EIS); Selectable Floating; Second Sense (S2); Analog Scan;

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