Application Note AN-RA-006

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

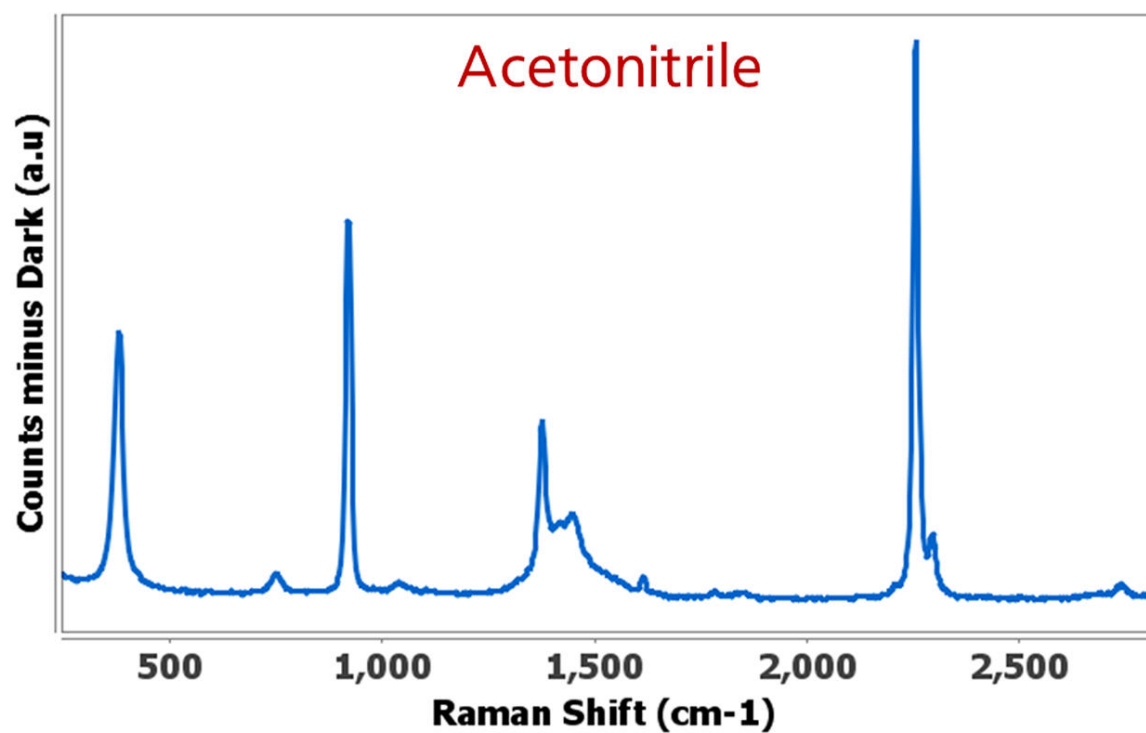
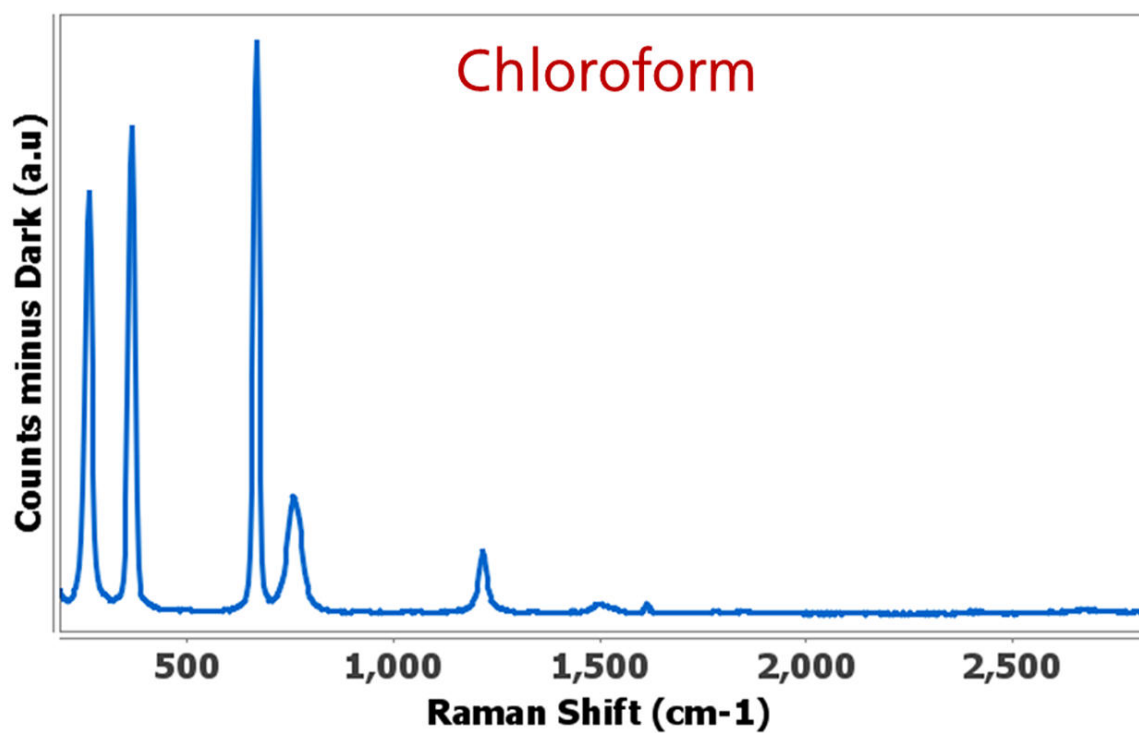
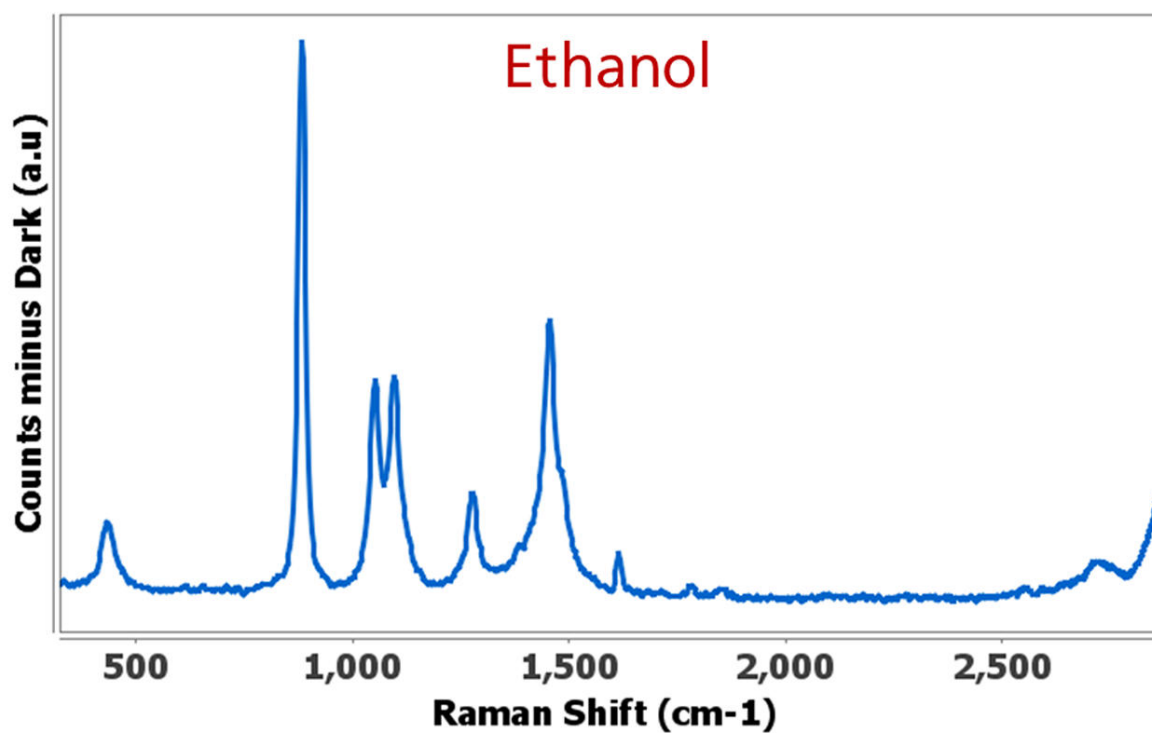
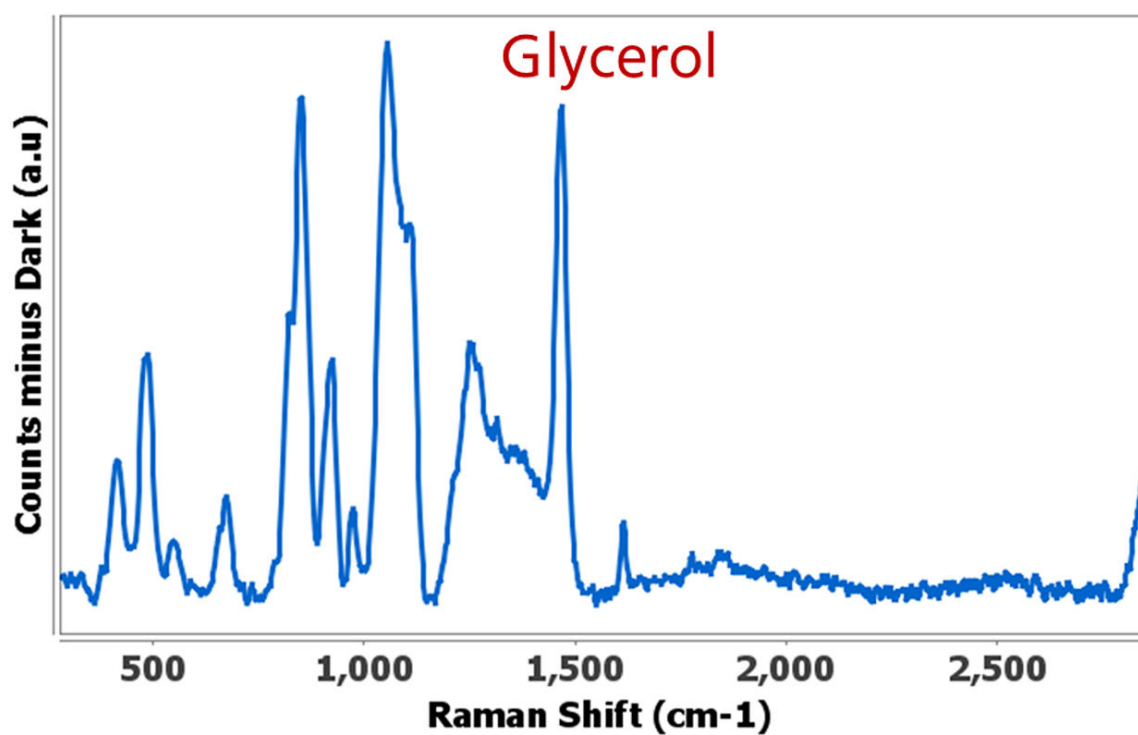
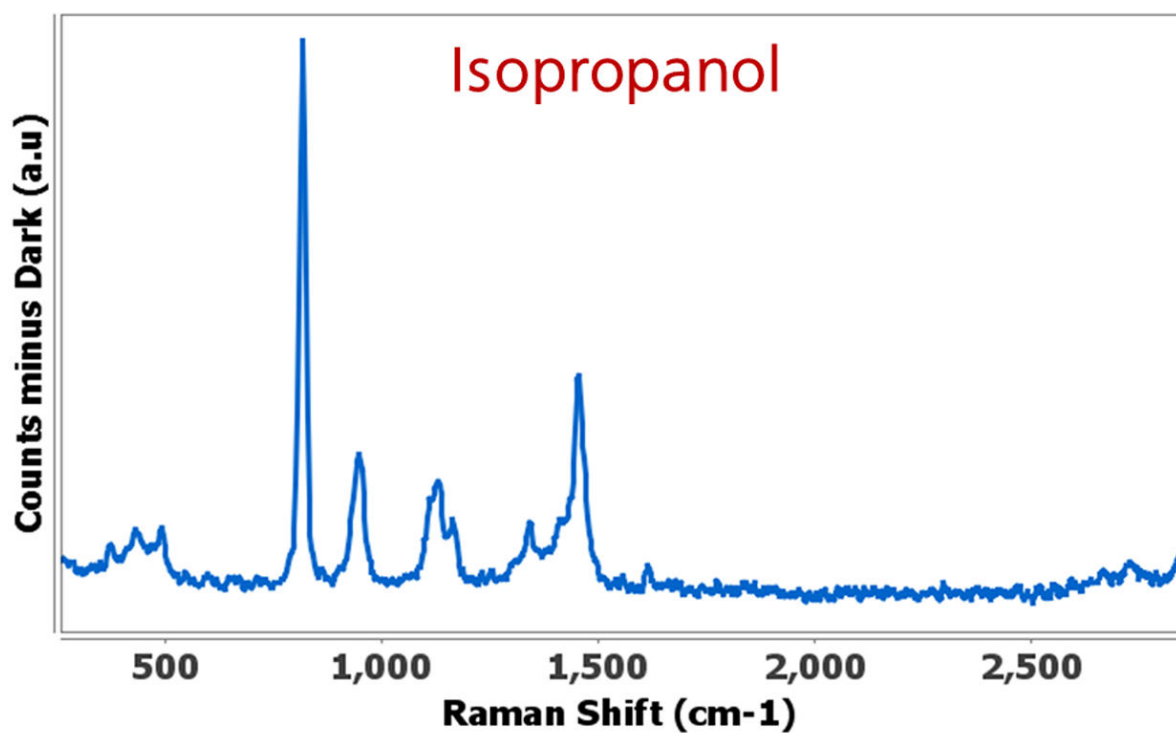
j)

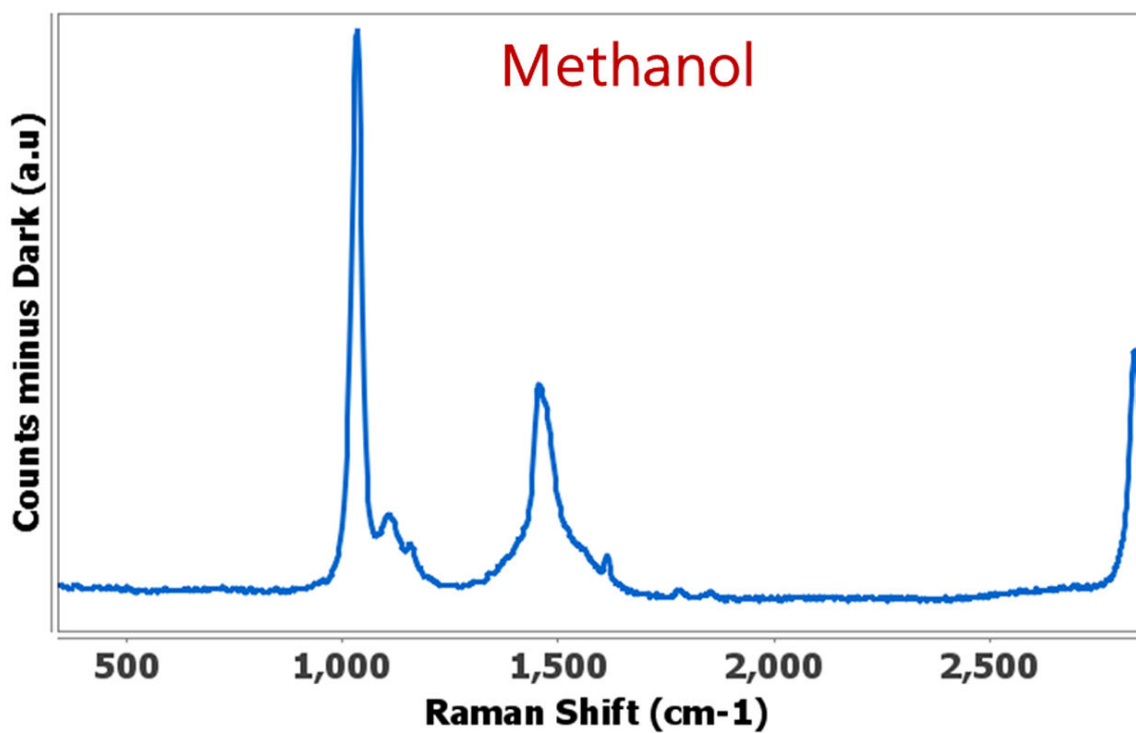k)

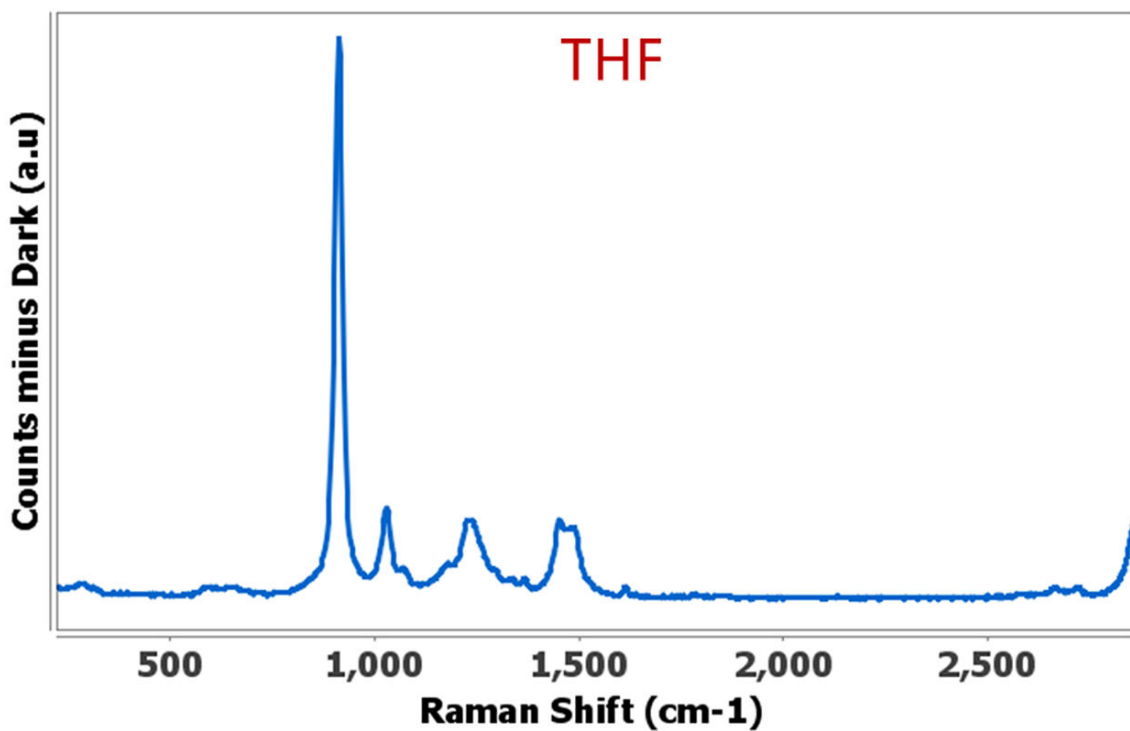l) Characteristic Raman spectra of (a) acetone, (b) acetonitrile, (c) cyclohexanone, (d) chloroform, (e) dichloromethane, (f) dimethylformamide (DMF), (g) dimethyl sulfoxide (DMSO), (h) ethanol, (i) glycerol, (j) isopropanol, (k) methanol, and (l) tetrahydrofuran (THF).

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
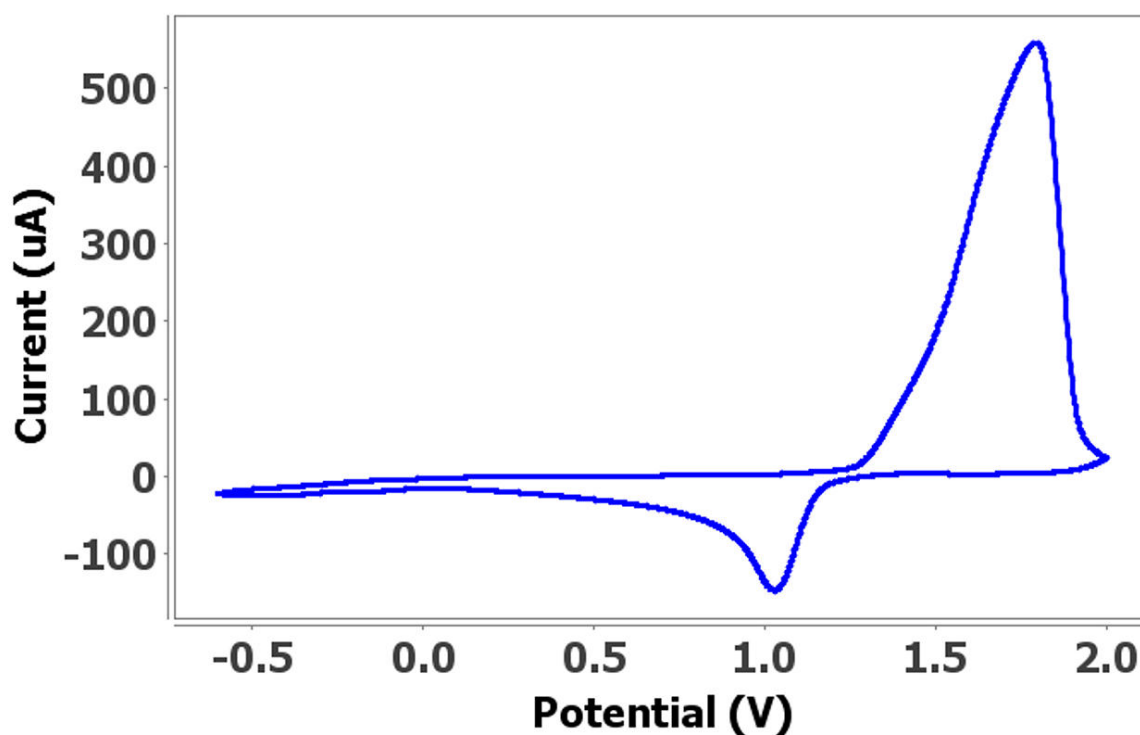
a)

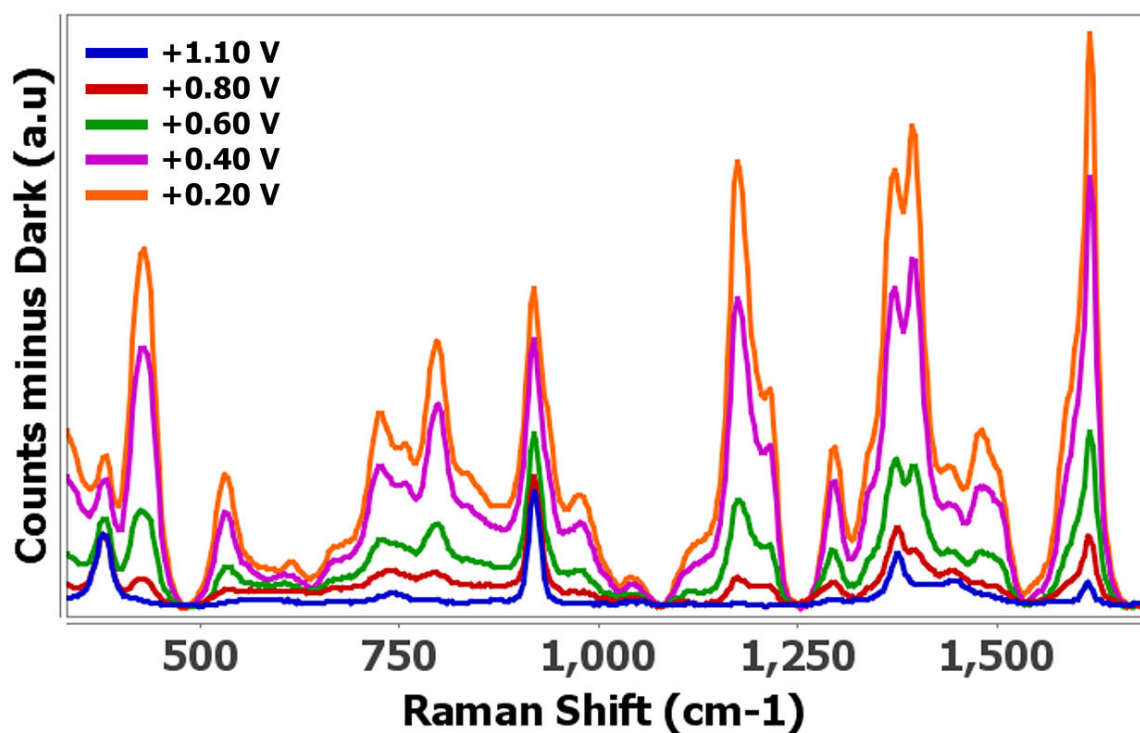b) a) Cyclic voltammogram and b) evolution of Raman spectra at different potentials obtained in 0.1 mmol/L crystal violet and 0.1 mol/L TBA in acetonitrile while scanning the potential from +0.70 V to +2.00 V and back to +0.60 V. The integration time was 2000 ms.

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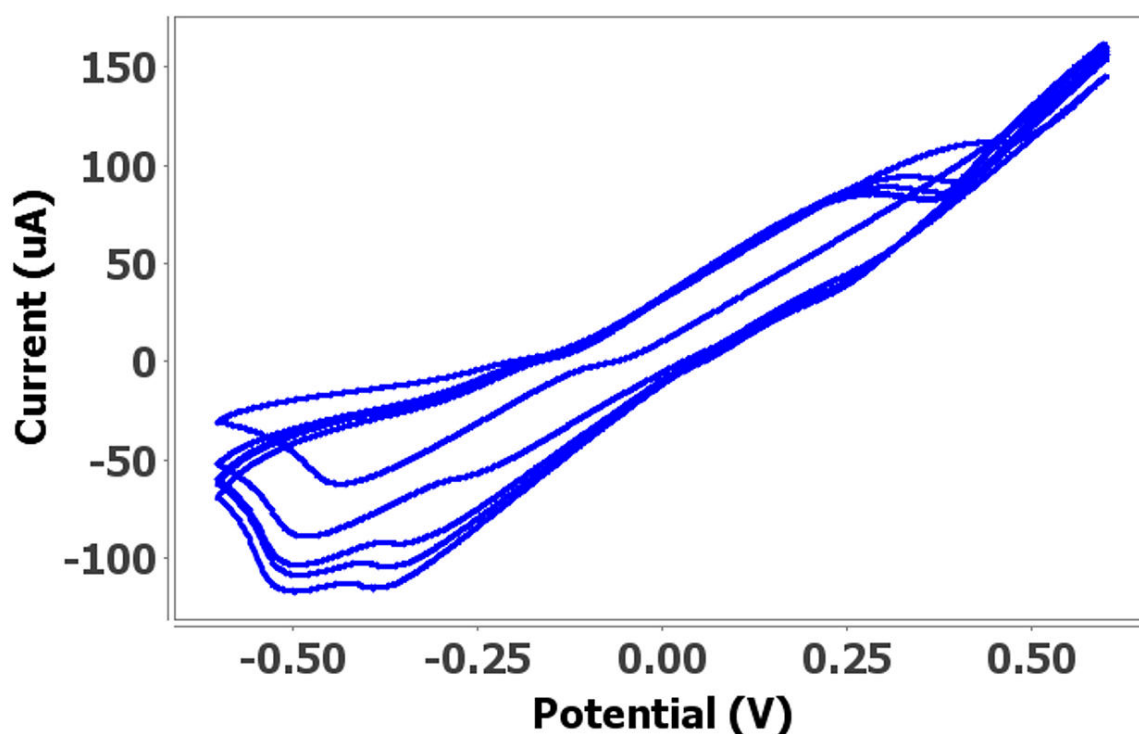
a)

b) a) Cyclic voltammogram and b) Raman spectra obtained in 0.5 mmol/L mancozeb and 0.1 mol/L TBA in DMSO while scanning the potential from +0.60 V to -0.60 V for five cycles. The integration time was 2000 ms.

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1. González-Hernández, J.; Ott, C. E.; Arcos-Martínez, M. J.; et al. Rapid Determination of the 'Legal Highs' 4-MMC and 4-MEC by Spectroelectrochemistry: Simultaneous Cyclic Voltammetry and In Situ Surface-Enhanced Raman Spectroscopy. *Sensors* **2022**, *22* (1), 295.
<https://doi.org/10.3390/s22010295>.
2. Ibáñez, D.; González-García, M. B.; Hernández-Santos, D.; Fanjul-Bolado, P. Detection of Dithiocarbamate, Chloronicotinyl and Organophosphate Pesticides by Electrochemical Activation of SERS Features of Screen-Printed Electrodes. *Spectrochimica Acta Part A: Molecular and Biomolecular Spectroscopy* **2021**, *248*, 119174.
<https://doi.org/10.1016/j.saa.2020.119174>.

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[AN-SEC-001 - Spectroelectrochemistry: an autovalidated analytical technique – Confirm results via two different routes in a single experiment](#)

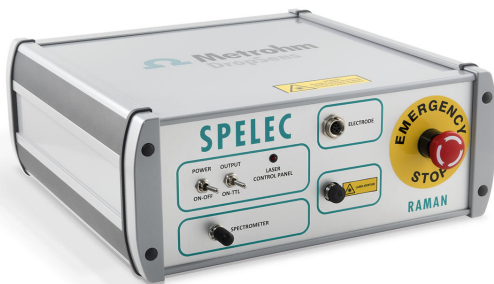
[AN-SEC-002 - Gathering information from spectroelectrochemical experiments – Calculation of electrochemical parameters from data](#)

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metrohm.jp@metrohm.jp

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
Spectroelectrochemical Raman instrument (785 nm laser)

SPELECRAMAN is an instrument for performing spectroelectrochemical Raman measurements. It combines in only one box a laser class 3B (785 nm $\pm$ 0.5), a bipotentiostat/galvanostat and a spectrometer (wavelength range 787 – 1027 nm and Raman shift 35 - 3000 cm⁻¹) and includes a dedicated spectroelectrochemical software that allows optical and electrochemical experiments synchronization.

Raman Probe

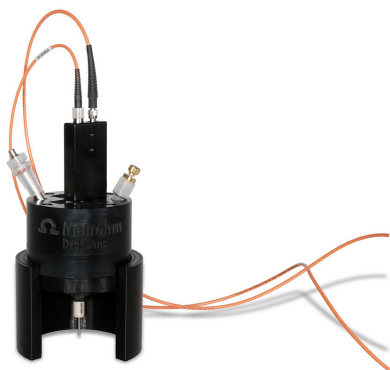Reflection probe designed to be used with a single excitation 785 nm wavelength (up to 500 mW). Suitable to work with DropSens Raman Cell for Screen-Printed Electrodes or with any conventional Raman set-up.

Raman spectroelectrochemical cell for conventional electrodes

Cell in PEEK for Raman spectroelectrochemical measurements. Designed be used with ref. RAMANPROBE and conventional Metrohm electrodes.