



Application Note AN-C-183

根据美国药典(USP)测定碳酸氢钾泡腾片口服混悬液中的钾

Method validation according to the U.S. Pharmacopoeia

Potassium bicarbonate effervescent tablets are used to prevent low levels of potassium in the blood [1]. Pharmaceutical manufacturers and laboratories are obliged to use the monographs from the United States Pharmacopeia (USP) and the National Formulary (NF) to test drugs and formulations.

The USP has embarked on a global initiative to modernize many of their existing monographs. As an alternative to atomic absorption spectroscopy (AAS), ion chromatography (IC) with non-suppressed conductivity detection has

been approved by the USP as a validated method to quantify the potassium content in potassium bicarbonate effervescent tablets for oral solution [2].

The Metrosep C 6 - 150/4.0 (L76) column provides the required separation of potassium. All acceptance criteria from the USP monograph «Potassium Bicarbonate Effervescent Tablets for Oral Solution» are fulfilled. The present IC method has been validated according to USP General Chapter <621> Chromatography [3].

SAMPLE AND SAMPLE PREPARATION

Potassium bicarbonate effervescent tablets for oral solution (Effer-K 25 mEq, potassium 978 mg, unflavored) from two different lots were evaluated in this application study.

Sample stock solutions with nominally 4890 mg/L potassium were prepared from at least 20 finely powdered tablets. A small amount (47.2 mg) of the resulting powder was transferred to a 2000 mL volumetric flask. 200 mL of ultrapure

water was added, and the flask was swirled until effervescence ceased. Then the solution was diluted to volume with ultrapure water and mixed well.

Sample solutions with nominally 15.0 mg/L potassium were prepared by transferring 1.533 mL of the sample stock solution to a 500 mL volumetric flask, then diluted to volume with ultrapure water and mixed well.

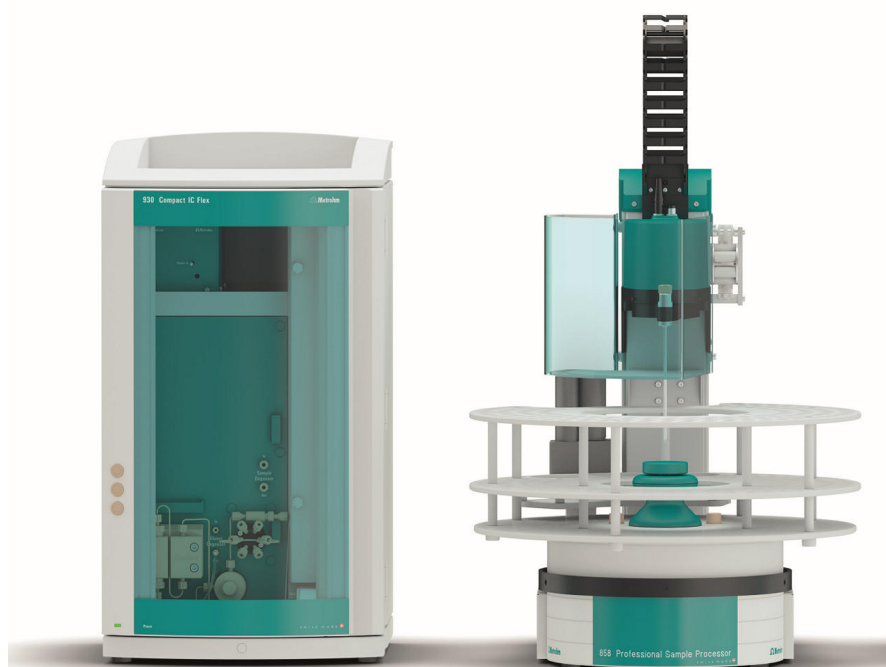


Figure 1. Instrumental setup including a 930 Compact IC Flex Oven/Deg and an 858 Professional Sample Processor.

EXPERIMENTAL

The working standard solution of 15 $\mu\text{g/mL}$ potassium was prepared from a certified 1000 $\mu\text{g/mL}$ potassium USP reference standard.

Samples and standard solutions were injected directly into the IC using an 858 Professional Sample Processor (**Figure 1**). Potassium was separated from all other cations using a

Metrosep C 6 - 150/4.0 column (L76) and the signal was recorded with a conductivity detector.

The calibration was performed using a single standard at 15 $\mu\text{g/mL}$ that was injected six times. The samples were analyzed in duplicate.

Table 1. IC method parameters as per USP monograph «Potassium Bicarbonate Effervescent Tablets for Oral Solution» [2].

Column with L76 packing	Metrosep C 6 - 150/4.0
Eluent	4 mmol/L nitric acid
Flow rate	0.9 mL/min
Temperature	30 ° C
Injection volume	20 µL
Detection	Direct conductivity

RESULTS

The IC assay for potassium content was validated according to USP Monograph «Potassium Bicarbonate Effervescent Tablets for Oral Solution» [2]. The accuracy of the potassium determination was calculated as 100% (Figure 2).

All acceptance criteria were fulfilled, e.g., asymmetry (tailing factors) for the potassium peak was <2, or the relative standard deviation (%RSD) of standard solutions was <0.5% (n=6) (Table 2).

Table 2. Required acceptance criteria as per USP monograph «Potassium Bicarbonate Effervescent Tablets for Oral Solution» [2].

Parameter	Actual	USP requirement	Status
% RSD standard solution (n=6)	0.05	NMT 0.5	Pass
Tailing factor	1.5	NMT 2.0	Pass
Resolution	3.89	NLT 3.0	Pass
Solution stability	0.08%	NMT 1.0%	Pass
Assay recovery	100.0%	90–110%	Pass
% RSD assay (n=6)	0.15%	NMT 1.0%	Pass

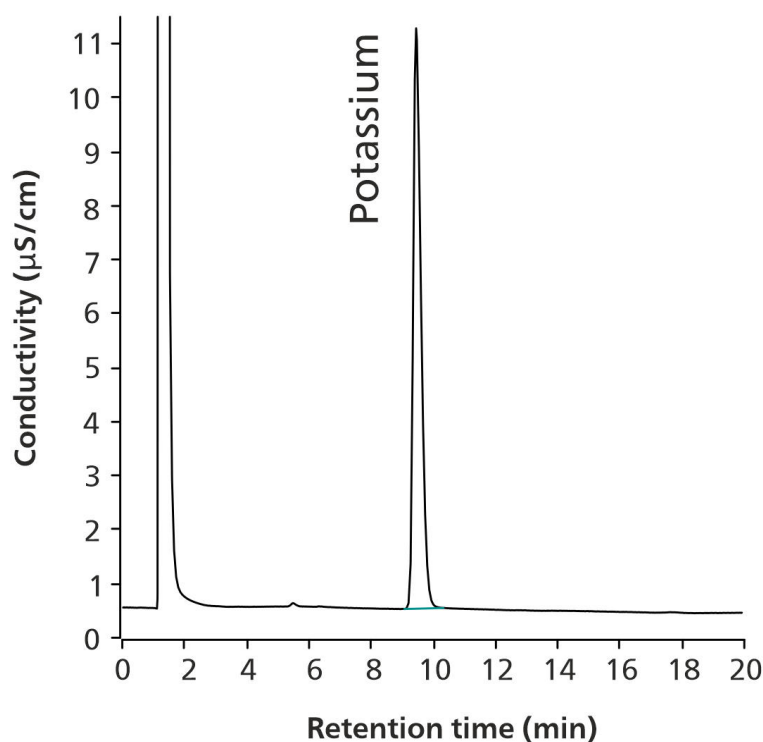


Figure 2. Chromatogram of 15.0 µg/mL potassium in sample solution (100.0% recovery of the nominal concentration).

CONCLUSION

The presented IC method for potassium in potassium bicarbonate effervescent tablets for oral solution with the Metrosep C 6 column (packing material L76) is officially included into the USP. Robustness and reliability of the

method was demonstrated following the guidelines of the USP General Chapter <621> [3]. The presented setup is suitable to quantify potassium according to USP requirements.

REFERENCES

1. Kardalas, E.; Paschou, S.S.; Anagnostis, P.; et al. Hypokalemia: a clinical update. *Endocrine Connections* **2018**, 7 (4), R135–R146.
<https://www.ncbi.nlm.nih.gov/pmc/articles/PMC5881435/>
2. U.S. Pharmacopeia. USP-NF Potassium Bicarbonate Effervescent Tablets for Oral Solution. *Monograph*.
https://doi.usp.org/USPNF/USPNF_M6719_4_02_01.html
3. <621> Chromatography.
https://doi.org/10.31003/USPNF_M99380_01_01

Internal references: AW IC IN6-1888-062018

CONTACT

瑞士万通中国
北京市海淀区上地东路1号
院1号楼7层702
100085 北京

marketing@metrohm.com.cn

CONFIGURATION



930 Compact IC Flex Oven/Deg

930 Compact IC Flex Oven/Deg 是智能型带柱加热炉、无抑制的 Compact IC 仪器,并且内置脱气装置。该仪器可使用各种分离和检测方法。


典型的应用范围:

- 阴离子和阳离子测定,无抑制的电导检测
- 使用 UV-VIS 或安培检测的简单应用



858 Professional Sample Processor – Pump

858 Professional Sample Processor – Pump 可处理体积在 500 μ L 至 500 mL 之间的样品。进行样品转移时,既可以使用内置的双向双通道蠕动泵、也可通过 800 Dosino 来进行。

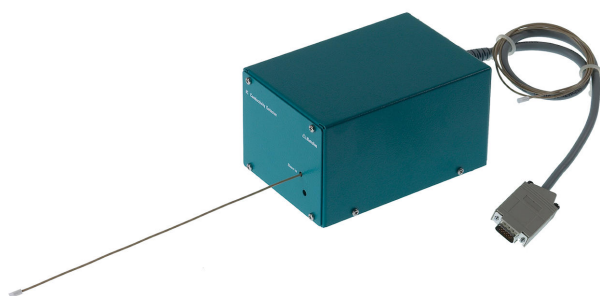

IC Conductivity Detector

用于智能型离子色谱仪的紧凑智能型高性能电导检测器。不凡的温度稳定性,受保护的检测器端子板内的整个信号处理过程以及新一代的 DSP(数字式信号处理)均能确保测量的准确性。归功于动态工作范围,无需进行范围变更(也不是自动进行)。