



Application Note AN-C-196

使用离子色谱法(IC)对三羟基甲基氨基甲烷(TRIS)进行纯度定量分析

Robust analysis with non-suppressed ion chromatography

Tris(hydroxymethyl)aminomethane (also known as TRIS, THAM, or tromethamine) is a common component of buffer solutions in the life sciences. It has a high buffering capacity between pH 7.2–9.0, a pK_a of 8.2 (20 ° C), and complexes with metal ions, making TRIS ideal for biochemistry and molecular biology applications [1]. TRIS buffers are used for DNA purification, separation of proteins with SDS-PAGE (sodium

dodecyl sulfate–polyacrylamide gel electrophoresis), or separation of nucleic acids with gel electrophoresis [2]. TRIS is also used to treat metabolic acidosis and can penetrate the cell membrane in its unionized form, therefore functioning as an intracellular buffer [3]. For these reasons, it is essential to control the purity of TRIS, especially for use in the pharmaceutical industry.

A robust isocratic ion chromatography (IC) method with a Metrosep C Supp 2 - 250/4.0 column and a methanesulfonic acid (MSA) eluent is ideally suited to determine TRIS and any cationic impurities. The microbore IC system (MB) is equipped with the IC Conductivity

Detector MB which is both sensitive and stable against MSA eluents. This guarantees low void volumes, long-term stability of the analytical system, and precise results for TRIS quantification.

SAMPLE AND SAMPLE PREPARATION

Samples were prepared from Trizma® base (TRIS) powder with p.a. quality (CAS 77-86-1, purchased from Sigma Aldrich No. 93350). For

method evaluation, two different concentrations of TRIS (10.37 mg/L and 103.7 mg/L) were dissolved in eluent (0.1% methanesulfonic acid).

EXPERIMENTAL

The microbore ion chromatograph 930 Compact IC Flex Oven/DEG/MB was equipped with the MSA-stable IC Conductivity Detector MB (**Figure 1**). An eluent consisting of 0.1% (v/v) MSA (15 mmol/L MSA) was used for this non-suppressed setup (**Table 1**). Samples were injected using the Metrohm intelligent Partial Loop Injection Technique (MiPT, **Figure 2**). This technique fills the 250 µL sample loop with a precisely measured and freely selectable volume (from 5 to 40 µL in this application study). During this process, a Dosino with a 2 mL Dosing Unit performs the precise dosing increments. MiPT enables calibration from a single standard, which was performed here in a range of 5140 mg/L TRIS.

The variable volume selection can also be applied to sample injection. In such situations, a small injection volume is selected, e.g., for a highly concentrated sample therefore omitting the manual dilution step.

Typical inorganic cations (i.e., lithium, sodium, potassium, magnesium, and calcium) were injected on the Metrosep C Supp 2 column to check for potential co-elution issues.



Figure 1. The IC Conductivity Detector MB shown here has a reduced cell volume and is inert against methanesulfonic acid.

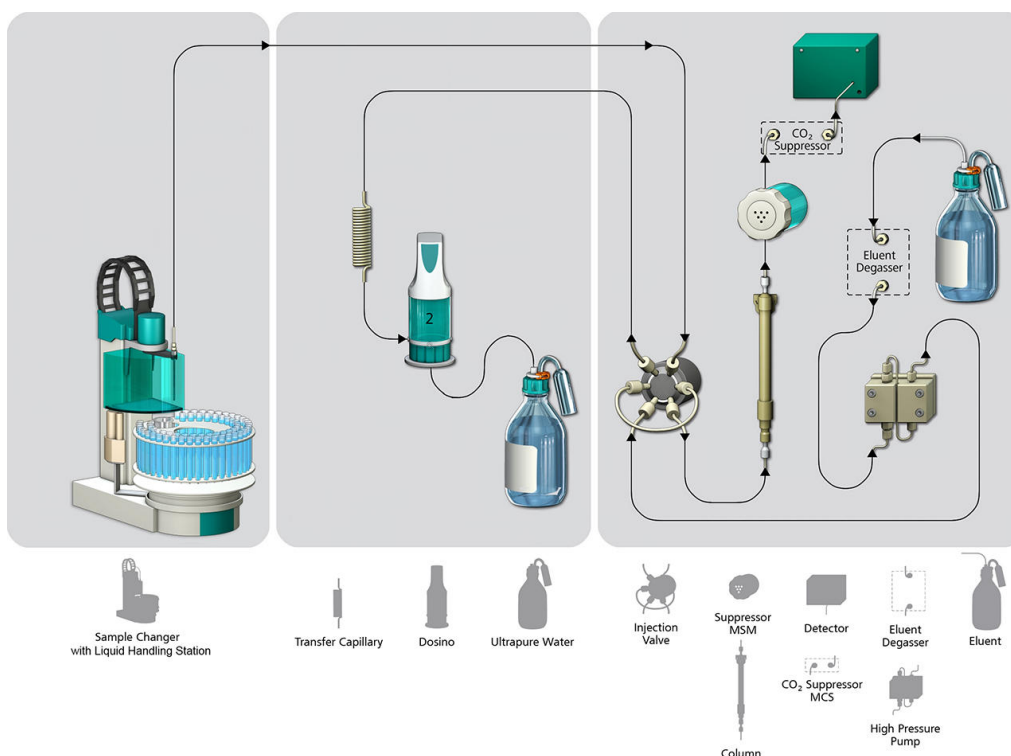


Figure 2. Illustration of the Metrohm intelligent Partial Loop Injection Technique (MiPT) flow path. With the help of the Dosino, the sample is transferred from the autosampler into a buffer loop to avoid contamination and carryover. The Dosino then precisely fills the sample loop with the desired injection volume in the μL range.

Table 1. IC method parameters for the microbore IC analysis of cationic impurities in TRIS.

Column	Metrosep C Supp 2 - 250/4.0
Eluent/diluent	$c(\text{MSA}) = 0.1 \text{ \% (v/v)}$
Flow rate	1.0 mL/min
Temperature	30 ° C
Injection volume	5–40 μL (MiPT)
Detection	Direct conductivity

RESULTS

TRIS determination is carried out in less than 8 min using isocratic elution on the MB IC system. The method was proven to be interference-free with regard to major cations as described above.

Sodium had a retention time of 4.1 minutes. Changing the method parameters, e.g., decreasing the column temperature to 20 ° C, will increase the resolution between sodium and TRIS. With the used method parameters (Table 1), precise determination is possible by using the peak height for evaluation. The recovery rates for 100 mg/L TRIS were 99–103% with a relative standard deviation of <3%, revealing the accuracy of this method.

CONCLUSION

Raw materials used in the pharmaceutical industry like solutions and buffers must fulfill the highest quality standards with respect to their exact concentration and purity.

The setup in this application study comprises a microbore IC system, an MSA-stable conductivity detector, and MiPT for automatic calibration

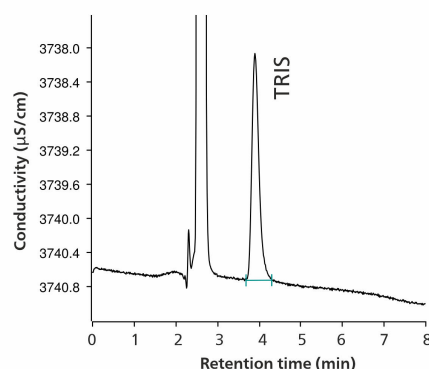


Figure 3. Chromatogram of 100 mg/L TRIS with 4 µL injection volume (MiPT).

with a single standard and flexible choice of sample injection volumes. The method is suitable for the quantification of TRIS in the range of 5200 mg/L. It guarantees robust determination of the common buffer component TRIS in an easy and precise way.

REFERENCES

1. Deutscher, M. P. *Guide to Protein Purification*; Gulf Professional Publishing, 1990.
2. Westermeier, R. *Electrophoresis in Practice: A Guide to Methods and Applications of DNA and Protein Separations*; John Wiley & Sons, 2016.
3. Sirieix, D.; Delayance, S.; Paris, M.; et al. Tris-Hydroxymethyl Aminomethane and Sodium Bicarbonate to Buffer Metabolic Acidosis in an Isolated Heart Model. *Am J Respir Crit Care Med* **1997**, 155 (3), 957–963.
<https://doi.org/10.1164/ajrccm.155.3.9117032>.

CONTACT

瑞士万通中国
北京市海淀区上地东路1号
院1号楼7层702
100085 北京

marketing@metrohm.com.cn

CONFIGURATION




930 Compact IC Flex Oven/SeS/PP/Deg/MB

930 Compact IC Flex Oven/SeS/PP/Deg 是智能型 Compact 离子色谱仪器,带有柱加热炉、序列抑制和用于抑制器再生的蠕动泵,以及内置的脱气装置。该仪器可使用任意分离和检测方法。

典型的应用范围:

- 阴离子或阳离子测定,序列抑制法及电导检测
- 针对微孔 (2mm) 应用经过优化,尤其适合耦合技术(IC-MS 或 IC-ICP/MS)

支持 MagIC Net 4.1 和以上版本

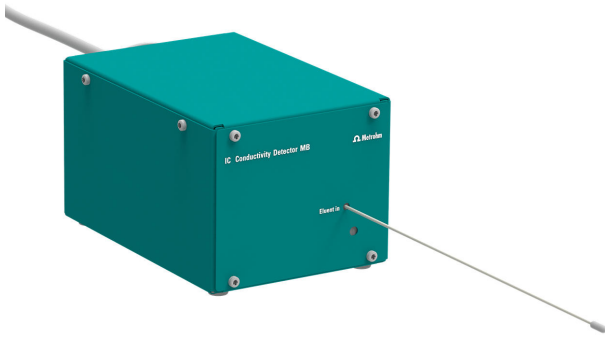

IC Conductivity Detector MB

用于智能型离子色谱仪的紧凑智能型高性能电导检测器。针对微孔柱经过优化。卓越的温度稳定性,在受保护的检测器端子板内完成整个信号处理过程以及最新一代的 DSP(数字式信号处理)均能确保测量的最高精确性。归功于动态工作范围,无需进行(也包括非自动)范围变更。

典型的应用范围:

- 阴离子或阳离子测定,化学抑制、序列抑制法或无抑制及电导检测
- 针对微孔 (2mm) 应用经过优化,尤其适合耦合技术(IC-MS 或 IC-ICP/MS)

规格说明书概览:

- 0 ... 15000 $\mu\text{S}/\text{cm}$ 无区段切换
- 测量池容量:0.3 μL
- 由不锈钢 X2CrNiMo17-12-2 (316 L) 制成的环形电极,与 MSA 兼容
- 最大工作压力:10.0 MPa (100 bar)
- 池温:20 ... 50 ° C,步幅为 5 ° C
- 温度稳定性: 0.001 ° C
- 基线噪音:° 0.2° nS/cm,是使用序列抑制法的典型值
- 毛细管:ID 0.18 mm

支持 MagIC Net 4.1 和以上版本



Metrosep C Supp 2 - 250/4.0

Metrosep C-Supp-2 系列长的分离柱是 Metrosep C Supp 2 - 250/4.0。Metrosep C-Supp-2 分离材料基于一种带羧基的聚苯乙烯-二乙烯苯共聚物。由于该分离材料经佳的钠/铵分离,故此该柱非常适用于除大量钠以外的小铵浓度测定。该柱可以搭配序列抑制法使用。故此,其特别适用于测定平均 $\mu\text{g}/\text{L}$ 范围及以下的浓度。