



Application Note AN-V-230

使用 scTRACE Gold 检测饮用水中的铬(VI)

Sensitive determination on the mercury film modified scTRACE Gold (DTPA method)

The guideline value for total chromium in the World Health Organization's (WHO) «Guidelines for Drinking-water Quality» is 50 µg/L. Chromium(VI) is more toxic than its trivalent form (Cr(III)) and is also less abundant. Therefore a robust and sensitive method is required to monitor its concentration in drinking water.

The mercury film modified scTRACE Gold can be used to monitor chromium(VI), offering easy handling and a high grade of stability. The sensor itself contains all necessary electrodes: the gold micro-wire working electrode, Ag/AgCl

reference, and the carbon auxiliary electrode integrated on a ceramic substrate which does not need intensive maintenance like polishing. Adsorptive stripping voltammetry (AdSV) using diethylenetriaminepentaacetic acid (DTPA) as a complexing agent and catalytic signal enhancement allows chromium(VI) detection at very low levels, with a limit of detection of 2 µg/L. The ability to replate the mercury film allows a quick and easy regeneration of the sensor. This method is best suited for field applications.

SAMPLE

Drinking water, mineral water, sea water

EXPERIMENTAL

Prior to the first determination, the ex-situ mercury film is deposited on the scTRACE Gold electrode. In the next step, the electrodes are cleaned with ultrapure water and the measuring vessel is emptied. The water sample, the supporting electrolyte with the complexing agent (DTPA) are pipetted into the measuring vessel. The determination of chromium(VI) is carried out with a 946 Portable VA Analyzer using the parameters specified in **Table 1**. The concentration is determined by two additions of a chromium(VI) standard addition solution. The scTRACE Gold is electrochemically activated prior to the first determination.

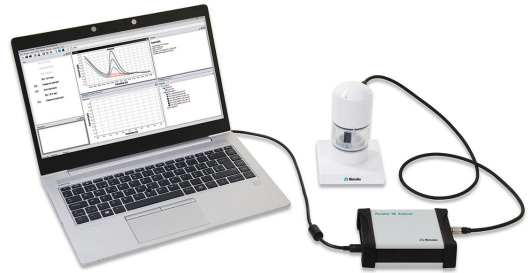


Figure 1. 946 Portable VA Analyzer

Table 1. Parameters

Parameter	Setting
Mode	DP – Differential Pulse
Start potential	-1.15 V
End potential	-1.65 V
Peak potential Cr(VI)	-1.4 V

ELECTRODES

- scTRACE Gold

RESULTS

The method is suitable for the determination of chromium(VI) concentrations up to 40 µg/L. The

limit of detection without deposition time is approximately 5 µg/L.

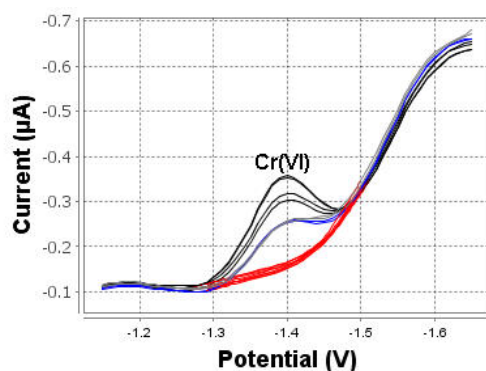


Figure 2. Determination of chromium(VI) in tap water spiked with 30 µg/L

Table 2. Result

Sample	Cr(VI) (µg/L)
Tap water spiked with 30 µg/L Cr(VI)	32.1

Internal reference: AW VA CH4-0596-042020

CONTACT

瑞士万通中国
北京市海淀区上地东路1号
院1号楼7层702
100085 北京

marketing@metrohm.com.cn

CONFIGURATION



946 Portable VA Analyzer (scTRACE Gold)

用于测定重金属,如痕量砷,汞,铜,铅,锌,镍,钴,铁、铋或锑之类重金属的便携式金属分析器。scTRACE Gold 适用的仪器版本。该系统由恒电位仪和集成了搅拌器与可更换式电极的独立测量台构成。该设备用 Portable VA Analyzer 软件驱动。电源由 USB 接口和内置的可充电电池提供。设备装在手提箱内交付,包含所有必需的附件。