



Application Note AN-CIC-034

通过燃烧离子色谱法(CIC)快速分析水中的可吸附有机卤素(AOX)

Measurement of AOCl, AOBr, AOI, and AOF according to DIN 38409-59 and ISO 18127

AOX (adsorbable organically bound halogens) is a complex parameter covering the sum of halogenated organic compounds adsorbable on activated carbon. Many of these organohalogenes and their degradation products pose serious risks to human health and the environment [1–4]. Monitoring them is essential to ensure appropriate water quality, to trace their sources, or to investigate the efficiency of AOX removal techniques in water treatment processes. Historically, AOX was determined via microcoulometric titration after adsorption of water samples on activated carbon and subsequent combustion (DIN EN ISO 9562 or

EPA 1650) [1,2]. By definition, based on the technical setup, AOX was comprised of adsorbable organically bound chlorine (AOCl), bromine (AOBr), and iodine (AOI)—but not fluorine (AOF)—as a sum parameter and not its individual fractions. Both **DIN 38409-59** and **ISO 18127** describe a validated procedure of adsorption and analysis via **combustion ion chromatography (CIC)** to determine AOCl, AOBr, AOI, AOF, and the sum parameter **CIC - AOX_(Cl)**. This Application Note explains the CIC method used to fulfill the requirements of these standards for AOCl, AOBr, AOI, AOF, and AOX analysis.

EXPERIMENTAL

This application is focused on the experimental approach of AOX analysis. More detailed information can be found in related Metrohm literature ([WP-081](#), [AN-CIC-033](#) – specifically about AOF). The complete validation dataset of DIN 38409-59 is available on the [Water Chemistry Society webpage](#).

The overall sample preparation procedure, i.e., preconcentration and adsorption of organically bound halogens, resembles that of DIN EN ISO 9562, as adsorption on activated carbon is a key point for both methods (**Figure 1**). While for AOF it is crucial that the samples are neutral to avoid adsorption of inorganic fluorine to the activated carbon, sample acidification is mandatory for the other organically bound halogens, similar to DIN

EN ISO 9562. For CIC-AOX_(Cl) determination (i.e., AOCl, AOBr, and AOI), samples need to be acidified with nitric acid to pH <2 prior to preconcentration (**Table 1**).

The adsorption of the organically bound halogens is handled in a semi-automated manner using the APU sim system from Analytik Jena (**Figure 1**). Two columns filled with activated carbon (at least 50 mg in each column) are connected in series and 100 mL of sample is passed through. The organically bound halogens adsorb to the activated carbon (using dedicated disposable columns for AOF and AOX determination, **Table 1**), while inorganic halogens are rinsed off (**Figure 1**).

After the semi-automated sample preparation is

finished, the complete content of the two adsorption columns is transferred into one or two separate ceramic boats for CIC analysis. Combustion occurs at temperatures above 950 °C in the presence of argon and oxygen (**Figure 1**). For pyrohydrolytic combustion, a water stream is essential as it converts the halogens into their hydrogenous forms. Chlorine, bromine, iodine, and fluorine are volatilized in the combustion step, transported into the absorber solution (ultrapure water) with an argon/oxygen gas stream, and transferred into the liquid phase (**Figure 1**). Dosinos guarantee precise automated liquid handling, e.g., the transfer of the aqueous sample into the IC for

analysis, or the water stream essential for pyrohydrolytic combustion.

The ion chromatographic separation is achieved on a Metrosep A Supp 5 - 250/4.0 column in combination with the A Supp 5 Guard/4.0. AOF (as F) elutes in under 7 minutes while AOX (i.e., Br, Cl, and I) elutes in less than 25 minutes (**Figure 2**). Automatic system calibration with MiPT (Metrohm intelligent Partial-Loop Injection Technique) is performed using inorganic anion standards for fluoride, chloride, bromide, and iodide (1 g/L standard solutions, TraceCert® from Sigma-Aldrich).

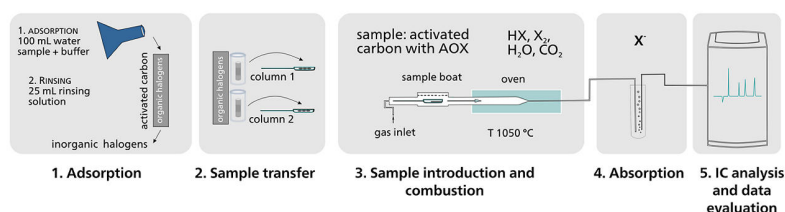


Figure 1. Schematic of the procedure for AOX (WP-081) and AOF (AN-CIC-033) analysis. The first step is adsorption performed with the APU sim (Analytik Jena) for semi-automated and standardized adsorption of up to six samples in parallel. After the second step of sample transfer into the combustion boats, the sample is automatically combusted (step 3, combustion module from Analytik Jena consisting of a combustion oven with Auto Boat Drive (ABD) and an autosampler (MMS 5000)). In the fourth step, the volatilized halogens are transported to the absorber solution via gas stream (920 Absorber Module). The last step (5) is the automatic analysis of AOB_r, AOCl, and AOI, or of AOF with the IC (930 Compact IC Flex) including data evaluation. The complete CIC process is fully automated and controlled by MagIC Net software from Metrohm.

Table 1. Parameters for AOF and AOX (AOCl, AOB_r, and AOI) sample preparation.

	AOF	AOCl, AOB _r , AOI
pH	Neutralized	Acidified to pH <2 with nitric acid
Buffer	0.5 mL 2 mol/L sodium nitrate	0.5 mL 2 mol/L sodium nitrate, acidified with nitric acid
Sample volume	100 mL	
	25 mL	
Rinsing solution	0.01 mol/L sodium nitrate	0.01 mol/L sodium nitrate, acidified with nitric acid
Absorption columns	Two activated carbon tubes (disposable, from Analytik Jena)	
	402-880.616	402-880.610
Flow rate APU sim	3 mL/min	

Performance checks of AOF and AOX determinations and the standard series for LOD determination (**Table 2**) are accomplished using organic reference standard solutions with varying concentrations (4-fluorobenzoic acid, 4-chlorobenzoic acid, 4-bromobenzoic acid, and 4-iodobenzoic acid), treated in the same way as

the samples.

As the procedure for the determination of AOX is rather complex, dedicated sample boats and charcoal (i.e., fluoride-free materials for AOF, **Table 1**) and blank measurements are essential to guarantee a low background and an appropriate blank correction (**Equation 1**).

RESULTS

Individual concentrations for AOCl, AOB_r, AOI, and AOF are calculated according to **Equation 1**. A sum parameter for AOX (CIC-AOX_(Cl)) is calculated using **Equation 2**. However, due to the novelty of this validated approach, CIC-AOX_(Cl) has not yet replaced AOX in water or wastewater regulations.

Dedicated materials and the sensitive analysis of the halogens with suppressed conductivity detection results in low blank values. Blank values were only measurable for fluoride and chloride (**Table 2**). The requirements for DIN 38409-59 are fulfilled—in fact, the overall procedure here is even more sensitive.

$$c(X_{ads}) = \left(c(X^-)_{IC} * \frac{V_{Abs}}{V_{Smpl}} \right) - \left(c(X_{BW}^-)_{IC} * \frac{V_{AbsBW}}{V_{SmplBW}} \right)$$

Equation 1.

$c(X_{ads})$	Mass concentration of individual adsorbable organically bound halogens (with X = Cl, Br, I, and F) in µg/L
$c(X^-)$	Halogen concentration in the sample's absorption solution in µg/L (with X = Cl, Br, I, and F) in µg/L
V_{Abs}	Final volume of the absorption solution in L
V_{Smpl}	Volume of the sample that was used for adsorption; always 0.1 L
$c(X_{BW}^-)$	Halogen concentration in the absorption solution of the blank in µg/L
V_{AbsBW}	Final volume of the absorption solution of the blank in L
V_{SmplBW}	Volume of the blank solution that was used for adsorption; always 0.1 L

$$c(CIC \cdot AOX_{(Cl)}) = c(AOCl) + c(AOBr) \cdot 0.4437 + c(AOI) \cdot 0.2794$$

Equation 2.

$c(CIC \cdot AOX_{(Cl)})$	Sum concentration of adsorbable organically bound halogens in µg/L as mass concentration based on chloride
---------------------------	--

During the DIN validation process, several water samples were analyzed from different

laboratories using similar setups (validation report: wasserchemische-gesellschaft.de).

Table 2. Blank, LOD (limit of detection), and DIN scope for the determination of adsorbable organically bound halogens. LODs are determined according to DIN 32645. For AOB_r and AOI, the LODs are determined using the calibration curve as no blank values were found. For AOF and AOCl, the blank method was applied (DIN 32645).

	Blank (µg/L)	LOD (DIN 32645) (µg/L)	Scope of DIN application (µg/L)
AOF	1.1	0.38	≥2
AOCl	2.6	1.36	≥10
AOBr	0	0.24	≥1
AOI	0	0.47	≥1

Using IC, it is now possible to not only determine the sum parameter $CIC-AOX_{(Cl)}$, but also to measure the individual fractions contributing to the AOX contents (**Figure 2**, [WP-081](#)) and to assess AOF ([AN-CIC-033](#)).

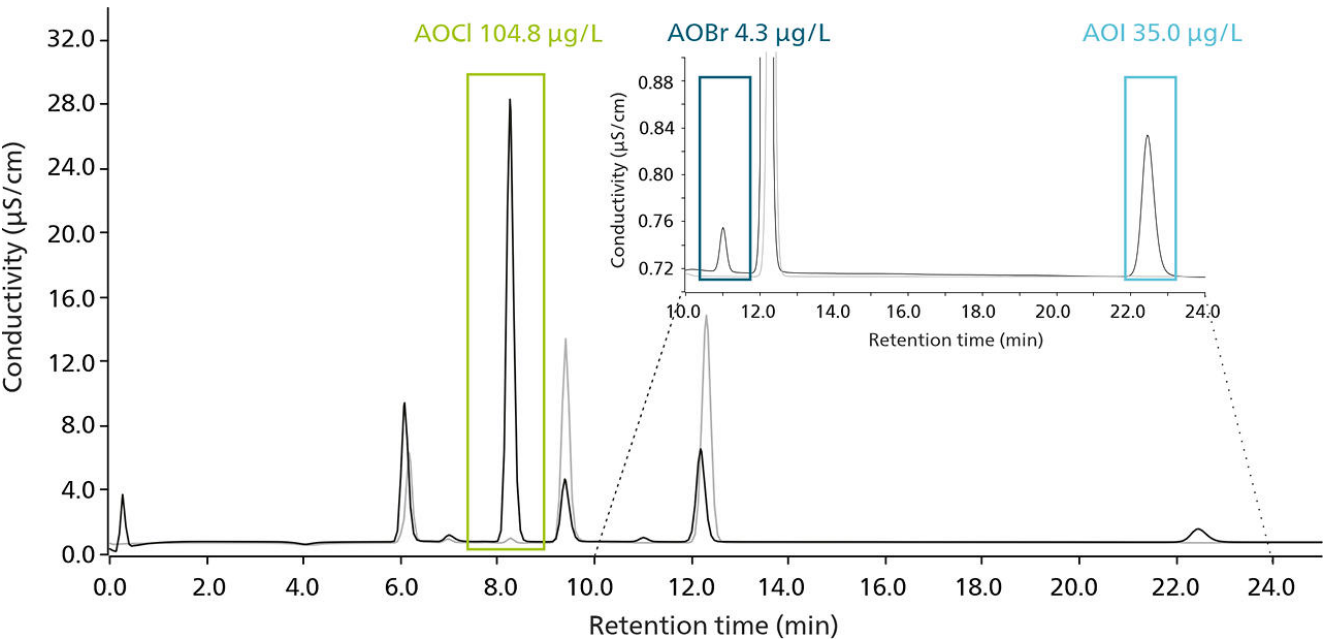


Figure 2. Chromatogram overlay of the blank and a wastewater sample for the determination of AOCl, AOB_r, and AOI measured from absorption column #1. To calculate the mass concentration of the individual AOX fractions, blank correction was performed according to Equation 1. No halogens were adsorbed on column #2, revealing the retention efficiency for AOX on column #1.

CONCLUSION

Overall, the validated procedure benefits from its easy, straightforward, and standardized handling, the precise determination of the analytes, the automatic calculation of results, and a low-maintenance, single-manufacturer setup.

A significant advantage of the standards DIN 38409-59 and ISO 18127 is that they allow the determination of adsorbable organically bound halogens as individual sum parameters, i.e., AOCl, AOBr, AOI, as well as AOF (a screening parameter for «total» PFASs, AN-CIC-033). Automation (e.g., automated eluent production, MiPT, intelligent and logical MagIC Net features)

improves repeatability, accuracy, and reliability of the results, saves valuable laboratory time for the liquid handling, standard, and eluent preparation, and allows 24/7 analysis – from which every laboratory, either research, routine, or governmental lab – can profit.

The world of organically bound halogens is so varied that these sum parameters enable insights about hot spots, transport pathways, but also particularly vulnerable regions in a very simple way, while complex targeted analysis, if at all, can resolve individual organically bound halogens for deeper investigations afterwards.

REFERENCES

1. Xu, R.; Xie, Y.; Tian, J.; et al. Adsorbable Organic Halogens in Contaminated Water Environment: A Review of Sources and Removal Technologies. *J Clean Prod* **2021**, 283.
2. Müller, G. Sense or No-Sense of the Sum Parameter for Water Soluble “Adsorbable Organic Halogens” (AOX) and “Absorbed Organic Halogens” (AOX-S18) for the Assessment of Organohalogens in Sludges and Sediments. *Chemosphere* **2003**, 52 (2), 371–379.
3. Dann, A. B.; Hontela, A. Triclosan: Environmental Exposure, Toxicity and Mechanisms of Action. *J Appl Toxicol* **2011**, 31 (4), 285–311.
4. Xie, Y.; Chen, L.; Liu, R. AOX Contamination Status and Genotoxicity of AOX-Bearing Pharmaceutical Wastewater. *J Environ Sci* **2017**, 52, 170–177.

Internal reference: AW IC CH6-1438-042021

CONTACT

瑞士万通中国
北京市海淀区上地东路1号
院1号楼7层702
100085 北京

marketing@metrohm.com.cn

CONFIGURATION



Metrosep A Supp 5 - 250/4.0

瑞士万通的分离柱以其高分离度可用于难度较大的分离任务。即便是复杂的分离问题也可用 Metrosep A Supp 5 - 250/4.0 轻松解决,且其解决方法可重现。该柱的高容量可在不进行样品前处理的情况下测定 150 mg/L 氯化物中的 1 µg/L 溴酸盐。此柱的使用范围远不止用来测定标准阴离子。在半导体工业中或发电厂的锅炉给水过程中控制高标准的纯度要求时,Metrosep A Supp 5 - 250/4.0 柱是适宜的选择。



Metrosep A Supp 5 Guard/4.0

Metrosep A Supp 5 Guard/4.0 对 IC 阴离子柱 Metrosep A Supp 5 和 7 进行可靠保护,使其不受样品或淋洗液的污染。

它含有与 Metrosep A Supp 5 相同的分离材料,也同样由 PEEK(聚醚醚酮)制成,并直接以几乎无死点容积的方式拧到相应分离柱上(«On Column Guard System»)。保护柱将延长分析用柱的使用寿命,且不会影响其色谱分离效率。建议使用 A Supp 5 Guard/4.0,其价格优惠且操作简单。



MiPT

用于安装 Dosino 以进行局部循环进样的附件组。



858 Professional Sample Processor – Pump

858 Professional Sample Processor – Pump 可处理体积在 500 µL 至 500 mL 之间的样品。进行样品转移时,既可以使用内置的双向双通道蠕动泵、也可通过 800 Dosino 来进行。



930 Combustion IC PP (AJ)

930 Combustion IC PP (AJ) 可使用在线燃烧分解(高温水解)方法分析各种可燃样品中的卤素和硫,然后测定离子色谱(Combustion IC)。它包括所有需要的部件,如 Analytik Jena 的 Combustion Module(2.136.0700)、920 Absorber Module、930 Compact IC Flex Oven/SeS/PP/Deg 和 MagIC Net 软件。必要时,930 Metrohm Combustion IC-Paket 可附加一台 Autosampler(Autosampler MMS 5000) 用于固体和液体样品。整套分析过程包括样品进入和样品分解均自动化进行,且由 MagIC Net 控制。



Autosampler MMS 5000 (AJ)

Analytik Jena(耶拿分析仪器公司)的 Autosampler MMS 5000 (AJ) 用于搭配 Metrohm Combustion IC 对液态和固态样品进行全自动分析。为了根据正确的样品类型对模块化的 Multi-Matrix 给样器进行调整,需使用液体用套件 (6.7303.000) 或固体用套件 (6.7302.000)。