



Application Note AN-BAT-003

用于锂离子电池的静电间歇滴定技术 (GITT)

The galvanostatic intermittent titration technique (GITT) is a powerful electrochemical method widely used for characterizing the kinetics and thermodynamics of battery materials.

GITT has three major applications for lithium-ion (Li-ion) batteries: the determination of the diffusion coefficient, open-circuit voltage (OCV) analysis, and overpotential/internal resistance analysis. Ultimately, this technique provides valuable insights into the electrochemical

behavior of materials under realistic operating conditions, making it an essential tool for researchers and engineers involved in battery development and optimization.

This Application Note presents an overview of GITT, detailing its principles, procedural steps, and applications in the field of energy storage. The following sections will outline the methodology and data analysis. INTELLO and its powerful suite of battery-oriented commands will be shown.

INTRODUCTION

A GITT procedure consists of a series of fixed current pulses that correspond to a known C-rate. Each current pulse is followed by a relaxation time, during which time no current passes through the cell. The current is negative during the discharge and positive during the

MEASUREMENT

Some assumptions are made during the analysis of data that influence the experimental conditions we apply to the cell during the measurement. The first is that each pulse step produces only a very small potential change. For that to be true, the current must be small relative to the capacity of the battery, which means that C-rates of $C/10$ and $C/20$ are common for the current pulses. The length of the current pulse is also kept relatively short, usually between 5 and 30 minutes.

The next assumption is that an equilibration state is reached during the relaxation step. Since the time for this to occur is different for every

cell and system, the length of the relaxation step varies more. In some cases it can be minutes, in others it can be longer—around 1–2 hours—and in extreme cases, more than 10 hours. It is important to take note of this when investigating new materials and to adjust the relaxation time accordingly. Since a complete GITT measurement requires that the battery is taken from fully charged to fully discharged and back again, it can take longer than a month in some cases to complete a full measurement. As a galvanostatic technique, it is the E vs. time signal which is analyzed at the end of the measurement.

In **Figure 1**, a typical GITT plot on a two-electrode battery is shown. There are two branches in the plot, corresponding first to when the cell is discharged to a lower cut-off voltage, and second to when the cell is charged to an upper cut-off voltage. An enlarged section of the discharge part of plot is shown in the insert of **Figure 1**. It's clear from this insert that the plot is made of a series of active (blue) and rest (orange) steps. Charging and discharging the battery in this manner, with a series of current pulses, has some interesting consequences that allow researchers to extract some extra information from the battery. These are described below.

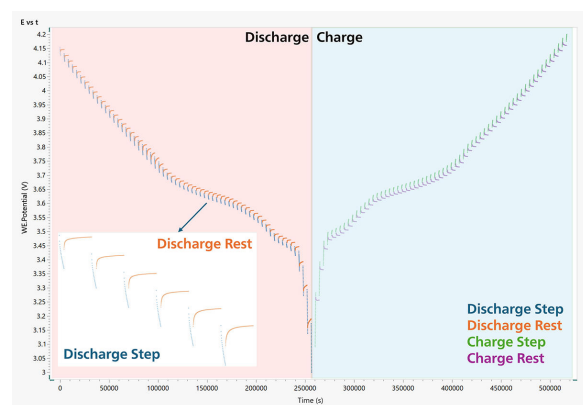


Figure 1. Complete GITT potential profile of a two-electrode battery being discharged and charged with current pulses that last 10 minutes, followed by a 60-minute rest period.

During a negative current pulse, the cell potential quickly decreases to a value proportional to the iR drop, where R is the sum of the uncompensated resistance R_u and the charge transfer resistance R_{CT} . Afterwards, the potential slowly decreases, due to the galvanostatic discharge pulse, in order to maintain a constant concentration gradient. When the current pulse is interrupted (i.e., during the relaxation time), the composition in the electrode tends to become homogeneous by Li-ion diffusion. Consequently, the potential first suddenly increases to a value proportional to the iR drop, and then it continues to slowly increase until the electrode is again in equilibrium (i.e., when $dE/dt \approx 0$) and the open circuit potential (OCP) is reached. Then, the galvanostatic pulse is applied again, followed by current interruption. This sequence of a discharge pulse followed by a relaxation time is repeated until the battery is fully discharged. During a positive current pulse, the opposite holds, which is repeated until the battery is fully charged.

Shown in **Figure 2** are the first five current pulses of the sequence; highlighted is the iR drop and ΔE_t – the voltage change during the current pulse, and ΔE_s – the change in the equilibrium or steady-state voltage due to the current pulse. All are useful in analysis of the GITT plot.

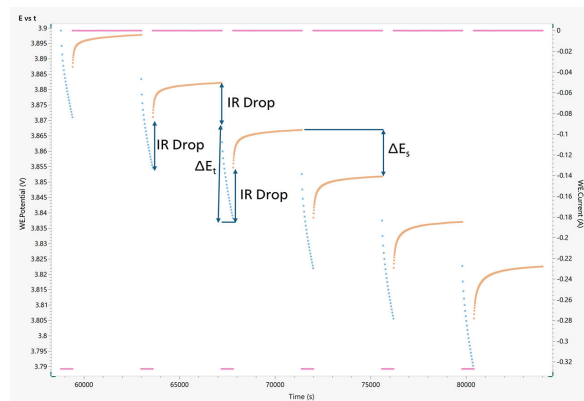


Figure 2. The first five current pulses are shown here. The pink bars across the top of the figure relate to the current applied to the cell during each step. Highlighted are the iR drop, the voltage change during the current pulse, and the voltage change in steady-state voltage due to the current pulse.

ANALYSIS

Diffusion Coefficients

The following equation is used to calculate the diffusion coefficient at each step (pulse plus relaxation) in the GITT procedure. Its derivation

from Fick's laws of diffusion is explained elsewhere [1–3].

$$D = \frac{4}{\pi} \left(\frac{iV_m}{z_A F S} \right)^2 \left[\frac{(dE/d\delta)}{(dE/d\sqrt{t})} \right]^2$$

Here, i (A) is the current, V_m (cm^3/mol) is the molar volume of the electrode, z_A is the charge number, F is Faraday's constant (96485C/mol), and S (cm^2) is the electrode area. Additionally, $dE/d\delta$ is the steady-state voltage change and $dE/d\sqrt{t}$ is the transient voltage change during

one galvanostatic titration step.

If sufficiently small currents (e.g., C/20) are applied for short time intervals (e.g., 10 minutes) then $dE/d\sqrt{t}$ can be considered linear. With these additional conditions, the previous equation can be simplified to:

$$D = \frac{4}{\pi\tau} \left(\frac{n_m V_m}{S} \right)^2 \left(\frac{\Delta E_s}{\Delta E_t} \right)^2$$

Here, τ (s) is the duration of the current pulse, n_m (mol) is the number of moles, V_m (cm^3/mol) is the molar volume of the electrode, S (cm^2) is the electrode area, ΔE_s (V) is the steady-state voltage change due to the current pulse, and ΔE_t (V) is the voltage change during the constant current pulse—eliminating the iR drop.

Typically, after the diffusion coefficient at each step is determined, it is then plotted as a

function of either the state of charge (SOC) or capacity of the battery. The changing state of charge is accompanied by physical changes in the electrode which can affect the diffusion of lithium ions. Monitoring the diffusion coefficient in this manner gives important insights into battery performance across the full charge/discharge cycle and helps researchers to optimize the performance of materials.

Open Circuit Potential

The OCP of a material at different states of charge is determined very accurately during the GITT procedure and contains some interesting thermodynamic information about the battery

material [1].

In this context, the OCP can also be defined as the chemical potential difference (μ) of lithium ions in the cathode and anode:

$$V(x) = \frac{\mu_{Li}^{Cathode}(x) - \mu_{Li}^{Anode}(x)}{e}$$

In this equation, x is the amount of lithium in the battery and e is the magnitude of the electronic charge.

Plotting this also as a function of SOC or capacity

is a useful tool to reveal changes in the electrochemical reaction of the battery as it is cycled.

Overpotential and Internal Resistance

In addition to the OCP, it can also be useful to look at the overpotential at each step. The overpotential is defined as the difference

between the measured cell voltage at the end of the current pulse (E_{meas}) and the voltage at the end of the relaxation step (E_{eq}) [1].

$$\eta = |E_{meas} - E_{eq}|$$

Considering the overpotential as a function of the SOC as well as the OCP can reveal kinetic and thermodynamic changes that are sometimes hidden by just looking at the overpotential in

isolation. It is also possible to instead consider changes in the internal resistance, which is essentially normalizing the overpotential to the applied current:

$$Internal\ Resistance\ (\Omega) = \frac{\eta}{i_{applied}}$$

Table 1. Details of the battery samples measured in this study.

Sample	Number of electrodes	Form	Capacity / mAh
1	Two	Cylinder	3270
2	Three	Pouch	4150

EXPERIMENTAL

To illustrate the principle of a GITT measurement, the technique was performed on two batteries with two and three electrodes, respectively, using VIONIC powered by INTELLO. There are two default procedures within the INTELLO library which are meant for GITT: one on a two-electrode battery and one on a three-electrode battery. The main parameters are C-rate, pulse, and rest durations, appropriate choices for which are discussed in the measurement section above. Also quickly

available are the voltage cut-offs, which are related to the battery under study (keep reading for some examples). For the 3-electrode procedure, cut-offs on both the WE-Second Sense (S2) and WE.potential signals are provided. Finally, parameters for adjusting the dynamic sampling rate are available. The standard plots are pre-built for convenience. Although we focus primarily on INTELLO in this Application Note, the same approach would be possible with an Autolab instrument and NOVA.

Table 2. Experimental details used in this Application Note.

Sample	C-rate	Pulse duration / min	Rest duration /min
1	C/10	10	60
2	C/10	10	10

In both cases studied, the samples were fully charged with a constant 0.5C before the GITT measurement was made. In the case of **Sample 1**, it was then discharged from 4.2 V to 3 V and

then charged again to 4.2 V. **Sample 2**, which has a different chemical makeup, was discharged from 3.65 V to 2.5 V and then charged again to 3.65 V.

RESULTS AND DISCUSSION

The GITT profile of **Sample 1** (the two-electrode battery) is presented in **Figure 3**. The conditions were satisfied such that it would be possible to calculate the OCP, overpotential, and internal resistance, and then plot this data vs. the SOC. However, since this is a two-electrode battery it is not possible to separate the diffusion contributions from the anode and cathode—therefore calculating the diffusion coefficient is impossible.

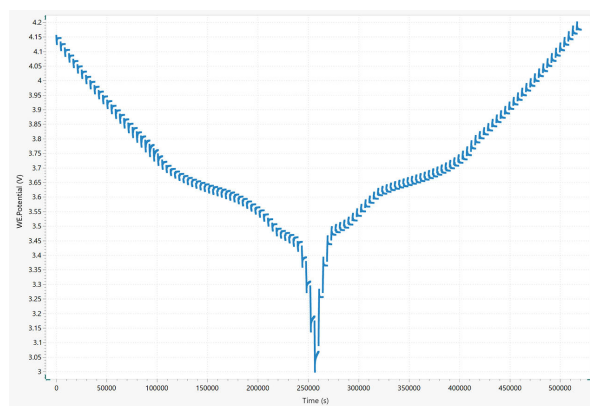


Figure 3. GITT profile of sample 1, a two-electrode battery with a capacity of 3270 mAh.

The GITT profile of **Sample 2** is shown in **Figure 4**. There are three potential signals in INTELLO: WE.potential, S2.potential, and WE-S2.potential. The standard connection to the three-electrode battery was made where the WE+S connected to the positive terminal, the CE+S2 connected to the negative terminal, and the RE connected to the reference electrode. The WE.potential (S to R) signal is the potential at the cathode, and the S2.Potential (S2 to R) signal is the potential at the anode. If the S2.potential is subtracted from the WE.potential, the voltage across the whole battery is calculated (i.e., the WE-S2 potential signal). Hence, in this case, the GITT profile is the WE-S2 potential vs. t. But the described analysis can also be performed independently on each potential signal, WE-RE and S2-RE, giving information about the diffusion of one or two half-cell materials.

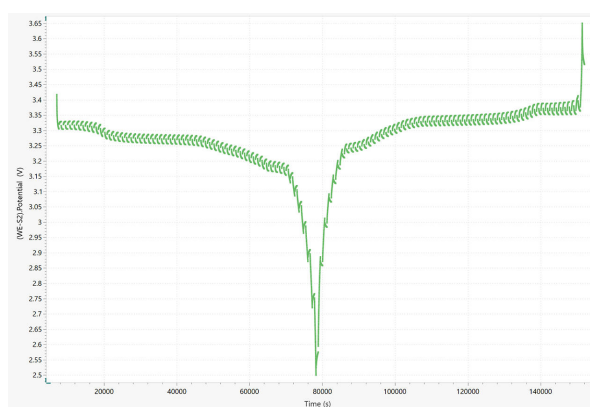


Figure 4. GITT profile of sample 2, a three-electrode battery with a capacity of around 4150 mAh.

Even though the experimental data would allow us to carry out further analysis and to calculate the diffusion coefficients, unfortunately it was not possible in this case due to the fact that the

surface area and the molar volume were not provided by the manufacturer of the battery used in this experiment.

REFERENCES

1. Kim, J.; Park, S.; Hwang, S.; et al. Principles and Applications of Galvanostatic Intermittent Titration Technique for Lithium-Ion Batteries. *J. Electrochem. Sci. Technol* **2021**, 13 (1), 19–31.
[DOI:10.33961/jecst.2021.00836](https://doi.org/10.33961/jecst.2021.00836)
2. Shen, Z.; Cao, L.; Rahn, C. D.; et al. Least Squares Galvanostatic Intermittent Titration Technique (LS-GITT) for Accurate Solid Phase Diffusivity Measurement. *J. Electrochem. Soc.* **2013**, 160 (10), A1842.
[DOI:10.1149/2.084310jes](https://doi.org/10.1149/2.084310jes)
3. Zhu, Y.; Wang, C. Galvanostatic Intermittent Titration Technique for Phase-Transformation Electrodes. *J. Phys. Chem. C* **2010**, 114 (6), 2830–2841.
[DOI:10.1021/jp9113333](https://doi.org/10.1021/jp9113333)

CONTACT

瑞士万通中国
北京市海淀区上地东路1号
院1号楼7层702
100085 北京

marketing@metrohm.com.cn

CONFIGURATION



Autolab PGSTAT302N

该恒电位仪/恒电流仪,具有 30 V 顺从电压,带宽为 1 MHz,可与我们的 FRA32M 模块联用,专门针对电化学阻抗谱而设计。

PGSTAT302N 是流行的 PGSTAT30 的后继款型。最大电流为 2 A,借助 BOOSTER20A 电流范围可扩展至 20 A,当电流范围为 10 nA 时电流分辨率为 30 fA。



Autolab DuoCoin Cell Holder

Autolab DuoCoin Cell Holder 具有 4 点开尔文镀金触头,可为您的电池研究提供较高精度的测量结果。一个通用附件可适应所有尺寸的标准电化学池以及较小和更大的非标准电化学池,可以一次处理二个电化学池。

Autolab DuoCoin Cell Holder 镀金触头和镀金印刷电路板可防止繁忙实验室中附件的腐蚀和损坏。

Autolab DuoCoin Cell Holder 具有可见电极标签和电缆连接件(与 Autolab 恒电位仪/恒电流仪电缆颜色相对应),可简化实验装置。Autolab 对细节的重视体现在 Autolab DuoCoin Cell Holder 底部的硅表面夹爪上(其可使复杂的实验装置具有稳定性)。



NOVA 是设计为通过 USB 接口控制所有 Autolab 仪器的软件包。

由电化学家针对电化学而设计,集成了超过二十余年的用户体验和最新的 .NET 软件技术,NOVA 使您的 Autolab 恒电位仪/恒电流仪拥有更强性能和灵活性。

NOVA 提供了以下的独特功能:

- 功能强大且灵活的程序编辑器
- 重要实时数据一目了然
- 强大的数据分析和绘图工具
- 集成化控制外围仪器,诸如万通 LQH 液体处理设备

[下载最新版本的 NOVA](#)



Autolab PGSTAT204

PGSTAT204 结合了小巧规格和模块化设计。该仪器包括基本恒电位仪/恒电流仪,其顺从电压为 20 V,最大电流为 400 mA 或 10 A,与 BOOSTER10A 组合使用。此恒电位仪可随时用附加模块进行扩展,例如 FRA32M 电化学阻抗谱(EIS)模块。

PGSTAT204 是一款经济实惠的仪器,可置于实验室的任何位置。具有模拟和数字输入/输出,可控制 Autolab 附件和外部设备。PGSTAT204 包括内置模拟积分器。与高性能的 NOVA 软件联用,可用于大多数标准电化学技术。