

区分不同制造商生产的异丙醇

This Application Note shows the rapid, non-destructive identification of isopropyl alcohol from two manufacturers using Raman spectroscopy following the creation of a suitable

library. The measurements with the handheld Raman spectrometer Mira M-1 require no sample preparation and provide immediate results that identify the samples unambiguously.

INTRODUCTION

Organic solvents like isopropyl alcohol are required in the production of many beauty and cosmetic products, paints, fragrances, and in synthesis reactions – the latter in particular in pharmaceutical applications.

In this work, isopropyl alcohol samples from two

different suppliers were analyzed and the differences in the spectra were investigated. Unexpectedly, the spectra showed some differences, allowing a supplier-specific identification.

EXPERIMENTAL

All spectra were measured using the Mira M-1 Raman spectrometer in auto-acquisition mode, i.e., integration times were determined automatically. A laser wavelength of 785 nm

and the Orbital-Raster-Scan (ORS) technique were used. The isopropyl alcohol samples were analyzed in vials using the vial holder attachment of the Mira M-1.

RESULTS AND DISCUSSION

Two different suppliers of isopropyl alcohol were evaluated. Due to the visible differences in the spectra indicating a possible contamination of

the samples the Raman spectroscopic distinction was possible (see **Figure 1**).

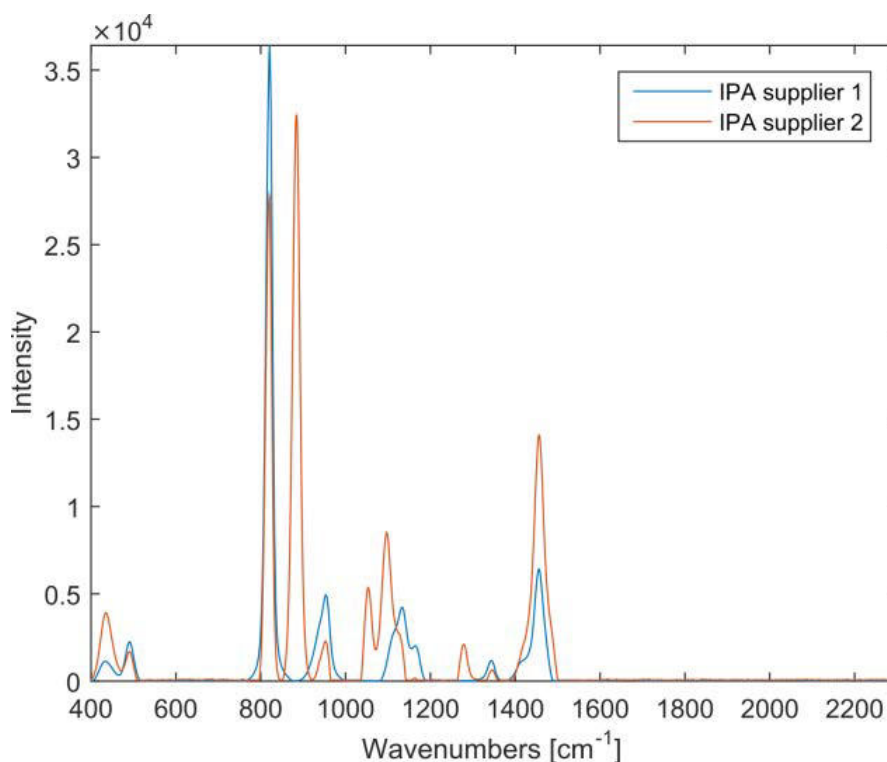


Figure 1. Spectra of isopropyl alcohol (Supplier 1 and 2) superimposed with a spectrum of ethyl alcohol

RESULTS AND DISCUSSION

The peaks at 1400–1470 cm^{-1} correspond to the asymmetric vibrations of $\delta(\text{CH}_2)$ and $\delta(\text{CH}_3)$. Between 600–1300 cm^{-1} , aliphatic chain vibrations for $\nu(\text{CC})$ can be found. In both samples of isopropyl alcohol, peaks in those areas can be observed.

Changes come up when focusing on the peaks at 883, 1049, 1095, and 1276 cm^{-1} , since those peaks, according to literature, suggest the presence of ethyl alcohol in the sample (see Figure 2).

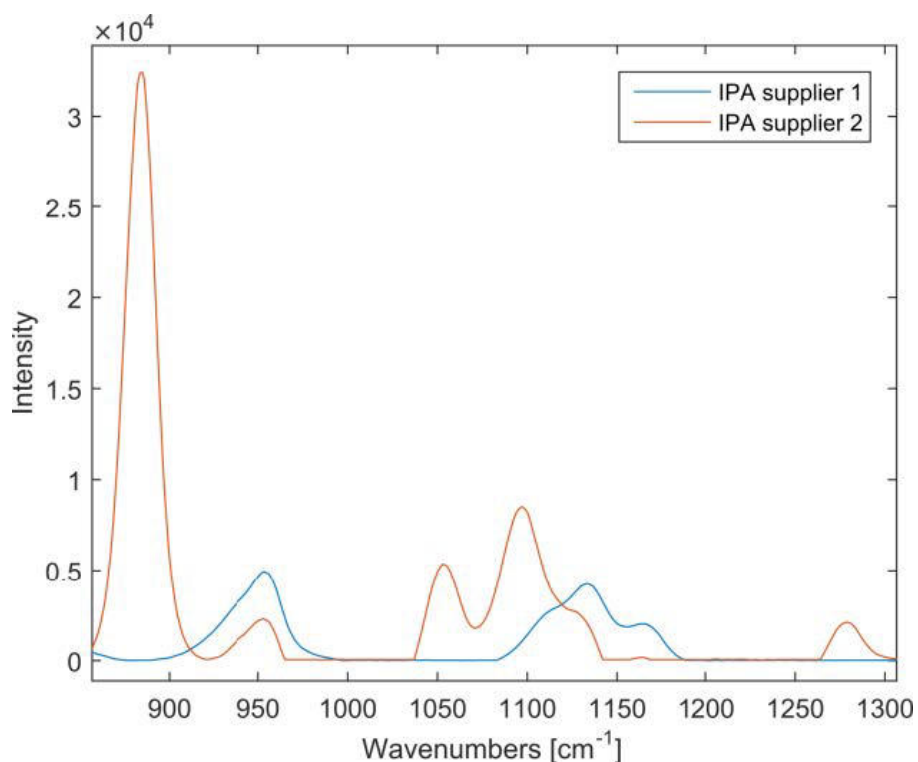


Figure 2. Differences between the isopropyl alcohol spectra.

CONCLUSIONS

Comparing the spectra of the two isopropyl alcohol samples, it becomes obvious that there is ethyl alcohol in the isopropyl alcohol sample of supplier 2. With Mira M-1, it was possible to

distinguish the different suppliers, proving that Mira M-1 is suited for the identification of incoming raw materials such as solvents and alcohols.

CONTACT

Metrohm Taiwan Ltd.
開封街二段 46 號 5 樓
108 台北市

info@metrohm.com.tw

CONFIGURATION



MIRA P Advanced

瑞士万通快速拉曼分析仪 (MIRA) P 是一款性能强大的手持式拉曼光谱仪,可用于各种材料的快速无损测定和验证,例如药物有效成分和赋形剂。MIRA P 小而坚固,配备了瑞士万通的ORS技术。MIRA P 符合 FDA 联邦法规 21 章第 11 款的规定

。Advanced Package 包含一个附加透镜,可用它直接分析材料或者在材料容器中分析(3b 级激光器),还有一个小管支架套筒用于分析玻璃小管中的样本(1 级激光器)。